



# Viewing Allergen Results

## Viewing Allergen Results on the EnSURE Touch

1. Open the Allergen App.

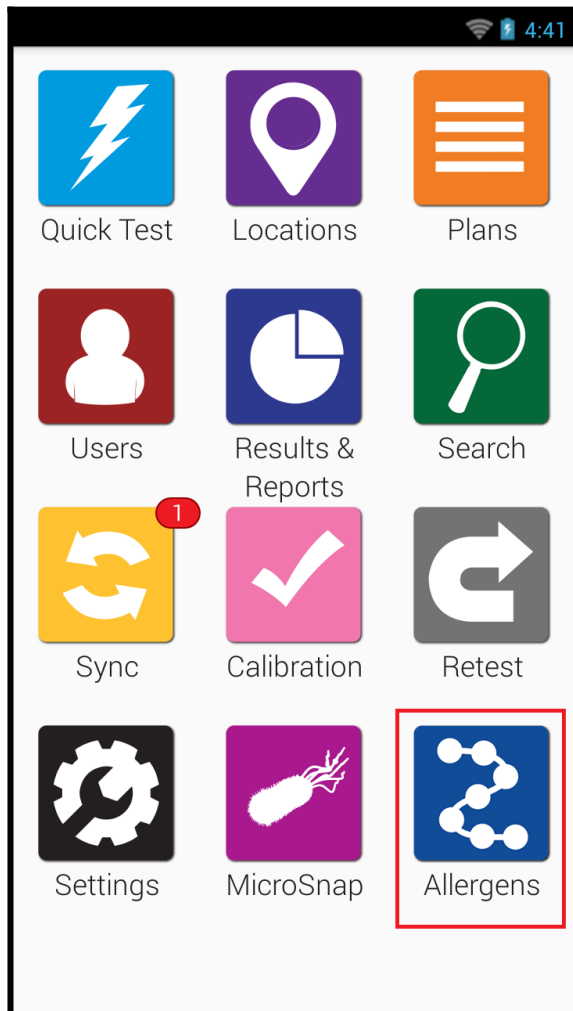

2. Tap Results.

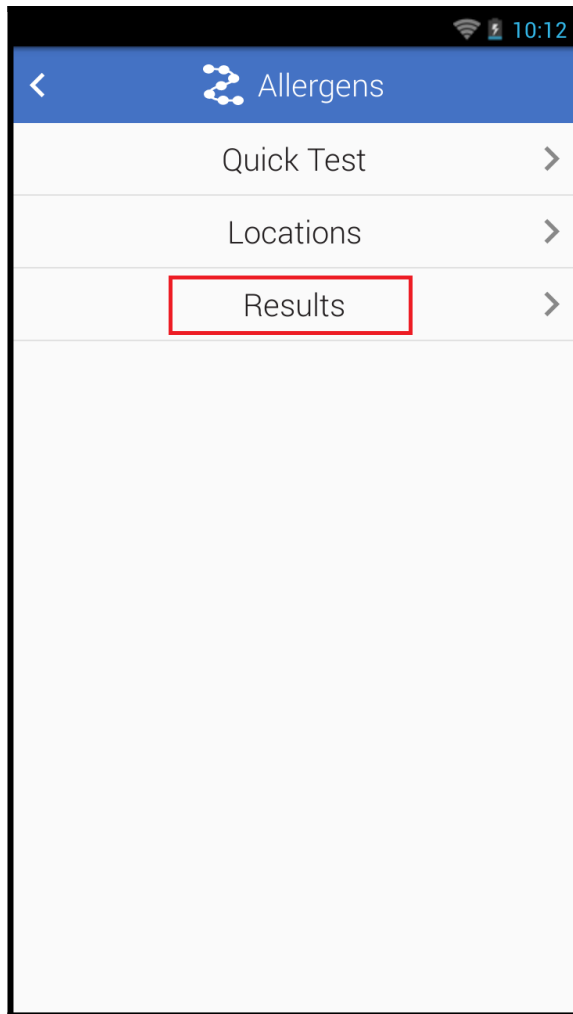

3. Tap To or From to a specific date range if required.

10:14

< Results ▾

From 10/20/2021

To 10/20/2021

Location

Results	%	Count	
✓ Pass	100%	2	>
⚠ Caution	0%	0	>
✗ Fail	0%	0	>
⚠ Invalid	0%	0	>

4. Tap the Location field to filter for specific locations.
5. Tap Pass, Caution, Fail, or Invalid to view a list of the results.

Results

From

10/20/2021

To

10/20/2021

Location

Results	%	Count	
✓ Pass	100%	2	>
⚠ Caution	0%	0	>
✗ Fail	0%	0	>
⚠ Invalid	0%	0	>