



Viewing Allergen Results

Viewing Allergen Results on the EnSURE Touch

1. Open the Allergen App.

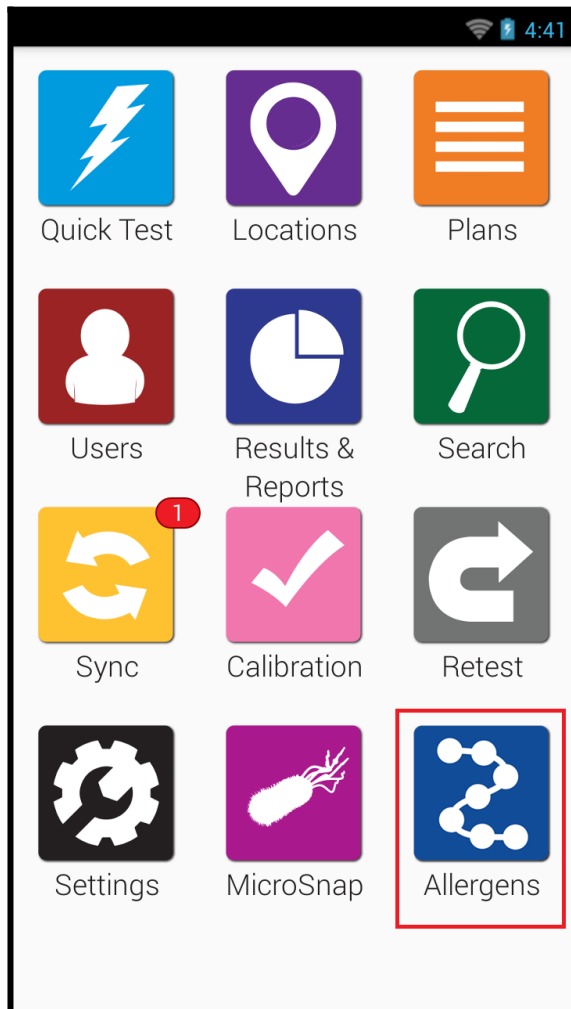

2. Tap Results.

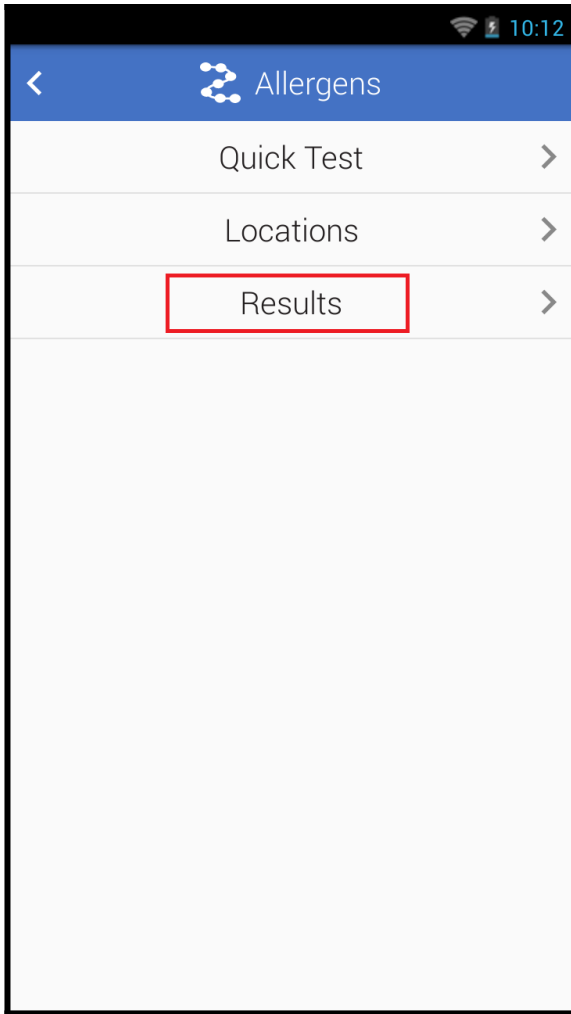

3. Tap To or From to a specific date range if required.

The screenshot shows a mobile application interface titled "Results". At the top, there is a blue header with a back arrow, a refresh icon, and the text "Results". Below the header, there are three filter fields: "From" (10/20/2021), "To" (10/20/2021), and "Location" (empty). These fields are enclosed in a red rectangular box. Below the filters is a table with the following data:

Results	%	Count	
✓ Pass	100%	2	>
⚠ Caution	0%	0	>
✗ Fail	0%	0	>
⚠ Invalid	0%	0	>

4. Tap the Location field to filter for specific locations.
5. Tap Pass, Caution, Fail, or Invalid to view a list of the results.



From

To

Location

Results	%	Count	
✓ Pass	100%	2	>
⚠ Caution	0%	0	>
✗ Fail	0%	0	>
⚠ Invalid	0%	0	>