

Download Center

In order to use this template effectively you may do one of two things to import your data into SureTrend:

1. Copy the headers from row 1 of this template to your exported SureTrend Excel file making sure you make any headers in your exported file match the ones from this template **exactly**. Spacing, punctuation and capitalization matter!
2. Remove the sample data from this template leaving only the headers in row 1. Copy and paste the data from your exported file into each column carefully. Be sure to be mindful of spacing, punctuation and capitalization as formatting matters! After inserting your data into this template you should be able to import it into SureTrend.

Using this template requires a minimum of SureTrend version 4.0.3.4 to use. It will not work with SureTrend 4.0.3.3 or earlier.

Need assistance?

Product support is included free of charge to *anyone!* Click the link below to open a new ticket and a specialist will be in touch with you soon.

You may also give us a call at [\(805\) 388-8007](tel:8053888007) and one of our sales or customer service associates can help you in no time.

[Open a ticket](#)

Download Details

Release date:	November 18, 2018
Last updated:	November 19, 2018
Current version:	1.0.0
Product type:	Excel Spreadsheet
File format:	.xlsx

File size:	21.30 KB
------------	----------

[Download Now](#)