

# Battery Replacement

## Introduction

This is a how to guide on switching out the battery on the **EnSURE Touch™**

1.) Disconnect the **EnSURE Touch** from the USB Power



2.) Remove the battery from the **EnSURE Touch**

To remove the battery, unscrew it with a small flat head screwdriver.



3.) Place the new battery in the **EnSURE Touch**



4.) Make sure it is flush to the unit when screwing the battery back in

5.) Let the **EnSURE Touch** fully charge (100%)



Your **EnSURE Touch** is ready to go!