


# Adding Zones to EnSURE Touch Locations

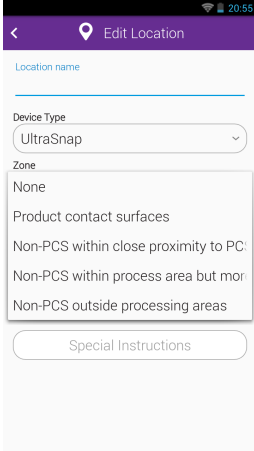

You can add zones to EnSURE Touch locations. This will allow data in the Results tables to be viewed by Zone type. The Zone is based on HACCP Zone criteria. There are 4 zones.

- Zone 1 = Product Contact Surface (PCS)
- Zone 2 = Non PCS within close proximity to PCS
- Zone 3 = Non PCS within process area but more removed from PCS
- Zone 4 = Non PCS outside processing areas

On the Ensure Touch, the Zone description is shown in the Zone text box. Zone detail can be added to locations either on the Sure Trend Cloud or on the EnSURE Touch. Before Zone detail can be added to locations the Zone feature needs to be enabled on the Sure Trend Cloud and Ensure Touch instrument.

## Enabling in SureTrend Cloud

1. Log into [SureTrend Cloud](#).
2. Click the **Sites** tab.
3. If necessary, click the site name to expand the site menu.
4. Click the **Site Settings** tab.
5. Scroll down and click **Location Settings** to expand the options.
6. Click the check box for **Enable Zone**.


Name	Logo	Active	
* Camarillo, US		<input checked="" type="checkbox"/>	

Locations	Plans	Instrument Users	SureTrend Users	Instruments	4 Site Settings
<b>Product Selections</b>					
<b>Security</b>					
<b>Plan Settings</b>					
<b>Location Settings</b> 5					
<input type="checkbox"/> Enable Group					
<input type="checkbox"/> Enable Surface					
<input checked="" type="checkbox"/> Enable Zone 6					
<input type="checkbox"/> Enable Line					
<input type="checkbox"/> Enable Cleaner					
<input type="checkbox"/> Enable Area					

## Enabling Zones on EnSURE Touch

1. From the **Home Screen**, tap **Settings**.
2. Tap **Customization**.
3. Tap **Locations**.

4. Tap the toggle to enable **Zones**.

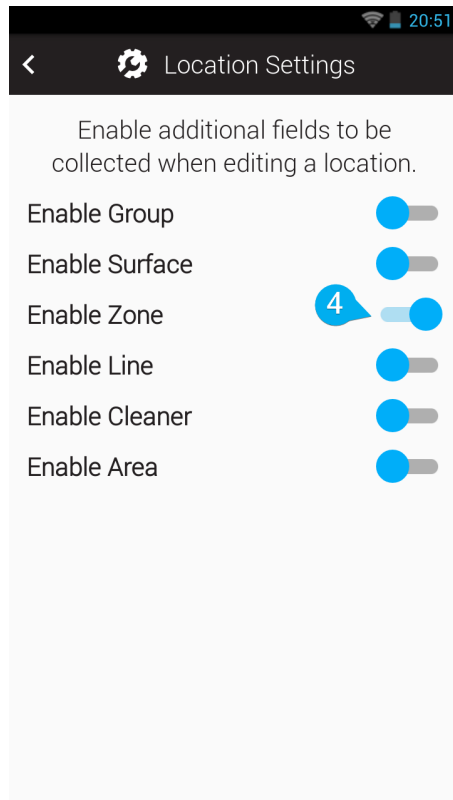

## Editing the Zone of Locations in SureTrend Cloud

1. Log into [SureTrend Cloud](#).
2. Click **Sites**.
3. If necessary, click your site name to expand the site menu.
4. On the **Locations** tab, click the **Edit** button for a location.
5. Enter a number 1-4 that represents the zone in the **Zone**.
6. Click **Save**.

Edit Location

×

Device/Assay/Test: \*

UltraSnap

▼

Location Name: \*

Hopper

Lower:

30

Upper:

60

Instructions:

Zone:

2

Group:

Select...▼

Surface:

Select...▼

Line:

Select...▼

Cleaner:

Select...▼

Area:

Select...▼

Image:

No Image

⬆

Global:

☒

6

Save

Cancel

## Editing the Zone of Locations on EnSURE Touch

1. From the **Home Screen**, tap **Settings**.
2. Tap **Locations**.
3. Tap the **Edit** button to enter the **Edit Locations**
4. Tap the **Edit** button for the location you want to edit.
5. Tap the **Zone** dropdown menu and select a zone option.

20:55

< Edit Location

Location name

Device Type

UltraSnap

Zone

None

Product contact surfaces

Non-PCS within close proximity to PCS

Non-PCS within process area but more

Non-PCS outside processing areas

Special Instructions

6. Tap the **Back Arrow** to return to the **Edit Locations**
7. Tap the **Back Arrow** again to exiting the **Edit Locations**.

**Contact Hygiena Technical Support for further assistance.**

- Phone: 1-888-HYGIENA (1-888-494-4362, option 2)
- Email: [techsupport@hygiena.com](mailto:techsupport@hygiena.com)
- [Submit a Support Ticket](#)
- [Schedule a Microsoft Teams meeting with support](#)