

Excluding Plans and Locations on EnSURE Touch

Each EnSURE Touch registered to the same site in a SureTrend Cloud account will share the same plans, locations, and users. If you want to have different plans and locations on each instrument while registered to the same site, you can use the Exclude feature to exclude specific plans and locations from appear on an EnSURE Touch.

Excluding Items from an Ensure Touch

1. Log into [SureTrend Cloud](#).
2. Click Settings and select [Instrument Management](#).
3. Click the serial number of your instrument.
4. Click the Locations, Plans, or Instrument Users tab.
5. Check the checkbox in the Exclude column to hide an item from the selected EnSURE Touch.

Instrument Management

Serial No	Instrument Name	Site	Calibration Check Status	Check Date	Pending Calibration	Model	Software	Last Sync
42967		Camarillo, US				EnsureTouch	1.140.9	9/22/2023, 4:11 PM

Instrument Serial No: 42967

Calibration Checks	Locations	Plans	Instrument Users	Power
	<div>Location Name</div> <div>Q</div> <div>ANES Cart</div> <div>ANES Keyboard</div> <div>ANES Machine</div> <div>ANES Mouse</div> <div>Arm Table</div> <div>B Scope Channel</div> <div>B Scope Distal</div> <div>Back Table</div>	<div>(All)</div>	<div>Exclude</div> <div><input checked="" type="checkbox"/></div> <div><input checked="" type="checkbox"/></div> <div><input checked="" type="checkbox"/></div> <div><input checked="" type="checkbox"/></div> <div><input type="checkbox"/></div> <div><input type="checkbox"/></div> <div><input type="checkbox"/></div> <div><input type="checkbox"/></div>	

Sync the Changes

1. On the EnSURE Touch, open the sync app.
2. Tap Sync now.
3. Wait for the sync to complete.
4. Check Plans, Locations, or Users to verify the items have been excluded from the instrument.

Contact Hygiena Technical Support for further assistance.

- Phone: 1-888-HYGIENA (1-888-494-4362, option 2)
- Email: techsupport@hygiena.com
- [Submit a Support Ticket](#)
- [Schedule a Microsoft Teams meeting with support](#)