
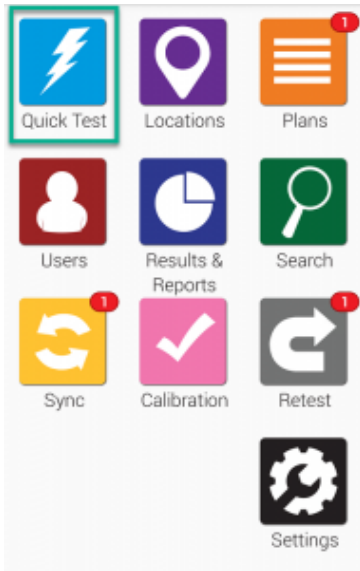


Performing a Quick Test

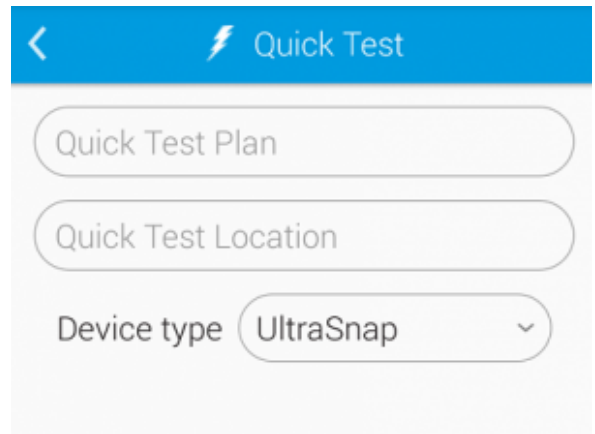

Click to enlarge

To get started, touch the **Quick Test** button from the Home screen. This will take you to the screen you'll see below.

The image shows a mobile application interface for a 'Quick Test'. At the top, there is a blue header bar with a white back arrow on the left and a white lightning bolt icon followed by the text 'Quick Test' on the right. Below the header, there are three rounded rectangular input fields. The first field is labeled 'Quick Test Plan'. The second field is labeled 'Quick Test Location'. The third field is labeled 'Device type' and has a dropdown menu with 'UltraSnap' selected and a small downward arrow on the right side.

[Click to enlarge](#)

Select a **Quick Test Plan** and a **Quick Test Location**. These will be from the collection of pre-defined plans and locations you've already created. If you do not have any plans or locations created yet you will be given a chance to create them on the Plans and Locations screens, respectively.

Select a **Device Type**.

If you have any custom fields for Tests active you may find more fields shown below Device Type such as Personnel, Room #, Inspection, and Product. You may enter your additional information into these fields before clicking **Run Test** to begin the Quick Test.