


# Turning Off Device Detection

## Introduction

Sometimes the device detection feature on EnSURE™ Touch will malfunction because the chamber in the instrument needs cleaning. Before you have time to thoroughly clean your instrument you can simply turn off the device detection if you have administrative rights.

## Step To Shut off Device Detection

1. On the home screen select settings

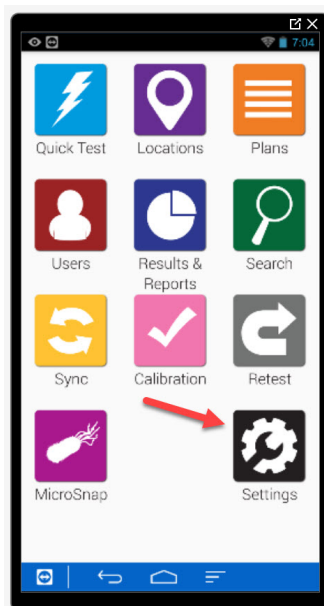


2. On the settings screen, select the General tab



3. On the General Settings tab select Auto-Detect



4. In the Auto-Detect Screen toggle the switch to off

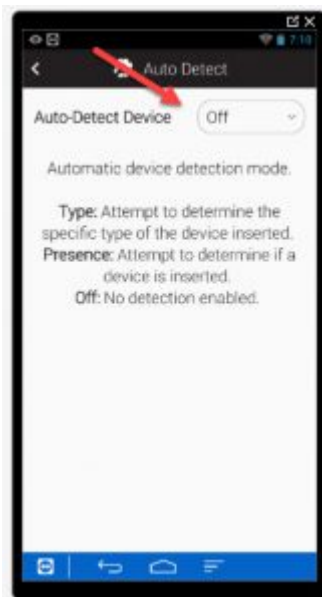

5. Return to the home screen.

You are now complete!

If you need further assistance please click on the icon for support to contact you.

