


Adding a Sample Location to SureTrend Cloud

Introduction

In this process, you will learn how to add locations to your SureTrend Cloud.

Steps

1. Log into your STC account

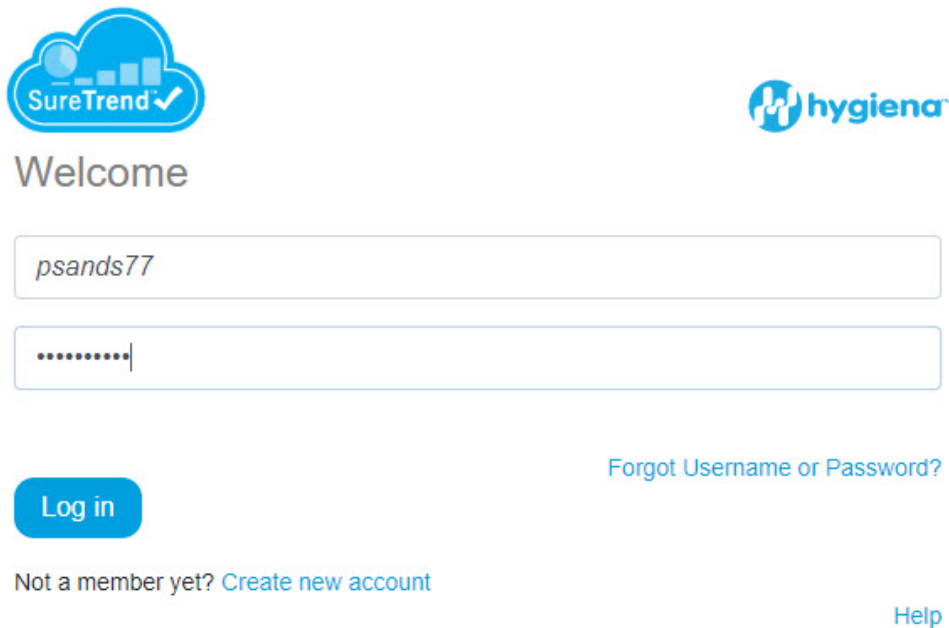

Welcome

psands77

.....

Log in

[Forgot Username or Password?](#)

Not a member yet? [Create new account](#)

[Help](#)









©2020 Hygienea, LLC

2. Click on the Instrument tab in the top left-hand corner




3. Select the instrument you wish to add the location to.

Instruments

	Serial No 	Instrument Name 
 	31576	Device 1
 	32253	Device 2
 	35624	Device 3

4. Click on the plus sign to add a new location.

Instrument Name: Device 2

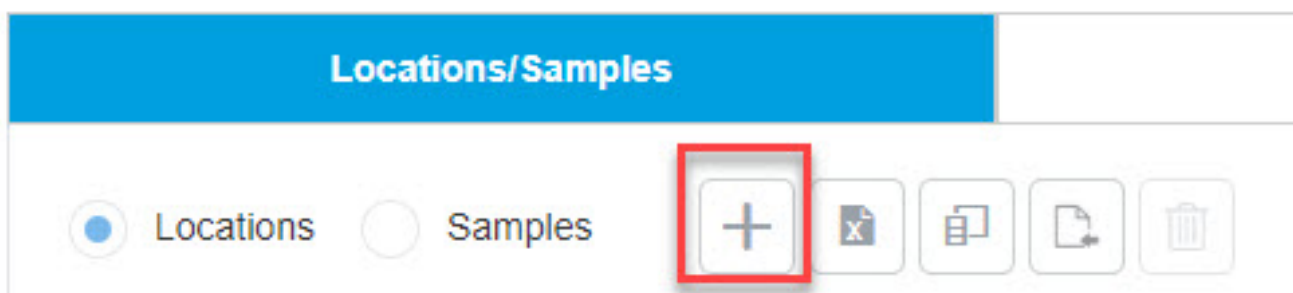

5. You can add the location that you need and adjust your thresholds. You can also add additional fields for further location detail. Some fields need to be enabled on the Ensure Touch.

New Location
✕

Location Name:

Lower:

Instructions:

Group:

Line:

Image:

No Image

Device Name:

Upper:

Zone:

Surface:

Cleaner:

7. Your new location is now added to all registered instruments. You can Exclude locations by clicking the exclude box. Locations can be edit by clicking on the pencil icon to make changes.

Locations
 Samples

		Location Name	Exclude	Image	Device Name	Lower	Upper
<input type="checkbox"/>	🔍		(All)		(All)	🔍	🔍
<input type="checkbox"/>		Kitchen	<input type="checkbox"/>		UltraSnap	65	120

Congratulations!

You are now complete! If you need further assistance please feel free to click on the link below.

