


Adding a New/Replacement EnSURE Touch to a Site


Introduction

This article will show you how to add a new or replacement EnSURE Touch to your SureTrend Cloud account.

Solution

1. Login to [SureTrend Cloud](#).
2. Navigate to the Instrument Tab.
3. Select  located on the right hand side above the Instruments table.
4. Enter an instrument name (optional), and select a Site. Note: If you only have a default site you will not need to select this.

If this is a replacement instrument be sure to select the proper site so your plans and locations will sync.

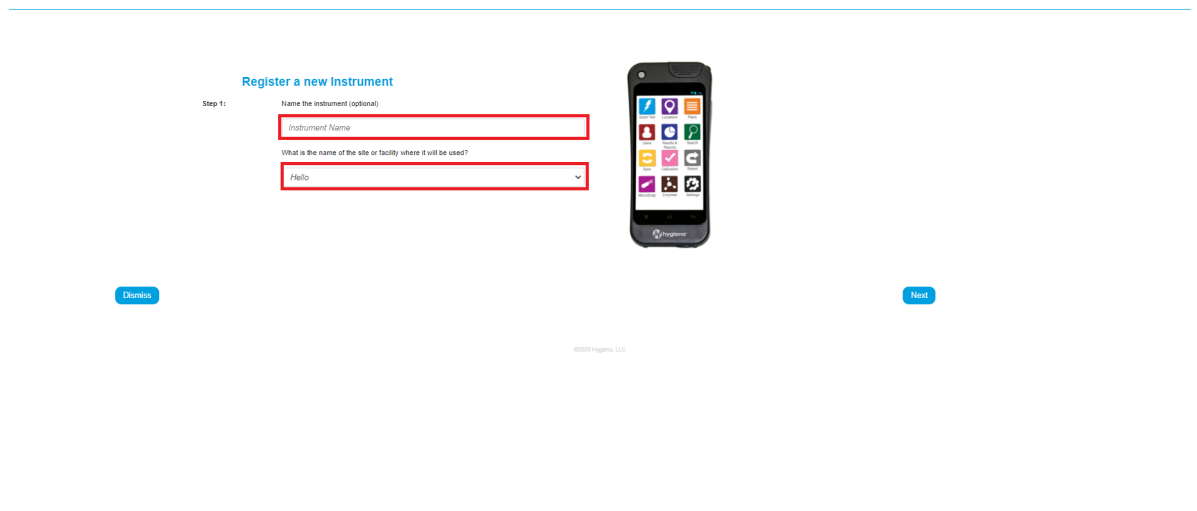


The screenshot shows a web interface for registering a new instrument. The title is "Register a new instrument". Below the title, it says "Step 1:". There are two input fields: "Instrument Name" (optional) and "What is the name of the site or facility where it will be used?". The second field has a dropdown menu with "Hello" selected. There are "Dismiss" and "Next" buttons at the bottom. A smartphone icon is shown to the right of the form.

5. On the EnSURE Touch tap the Sync App and when prompted enter the Username and Code.

Register a new Instrument

- Step 2: Connect Instrument to the Internet via Wi-Fi
- Step 3: On the Instrument, select the Sync icon on the home screen. Enter the Username and code on instrument and press the register button. Wait for status to change to successful
- Username:
- Code:
- Status: Pending

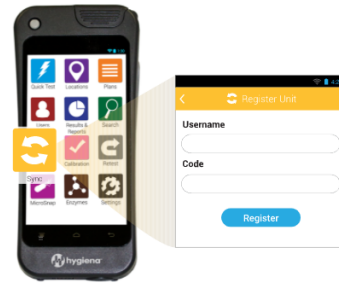

Cancel Registration

Done

6. Tap the sync now button to add your existing plans and locations.
Note: Plans and Locations are linked by the site the instrument are assigned to.
7. Your EnSURE Touch is now registered.