

Product Selection instructions for EnSURE Touch


Introduction

Product Selections allows you to choose what apps and test devices are available on the EnSURE Touch™. In addition to testing ATP the EnSURE Touch also allows you test Micosnap and Enzymes. This article will show you how enable or disable products from EnSURE Touch and SureTrend Cloud.


Solution

EnSURE Touch


1. Touch the **Settings** App from the home screen. Next, touch **Customization**.

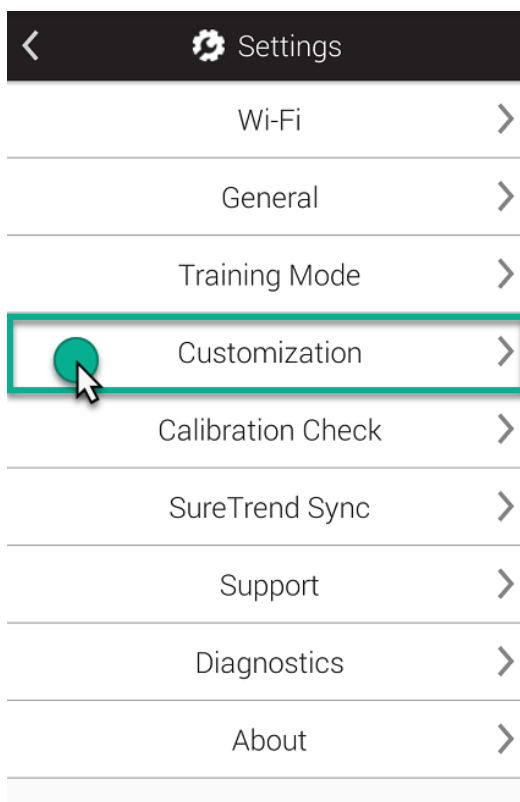

2. Select **Product Selections** □.

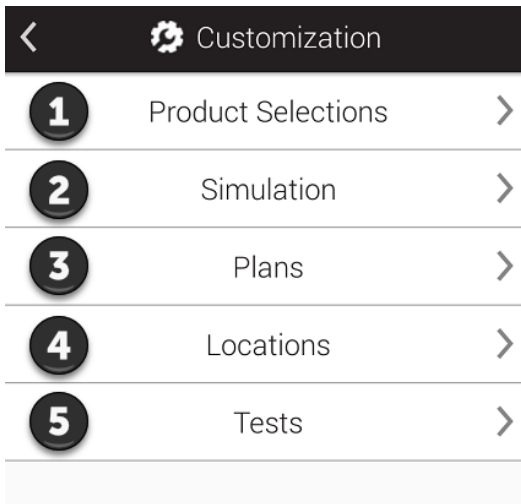

3. Select the product(s) you would like to enable/disable.

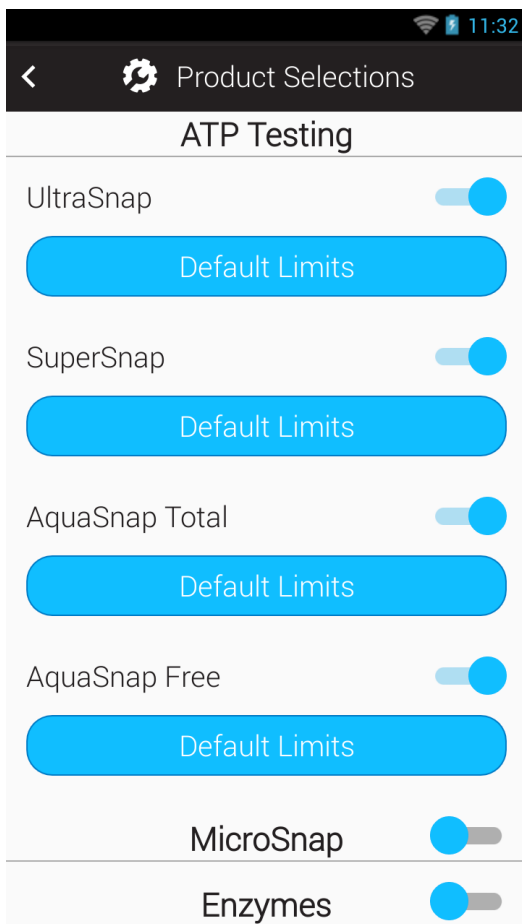

ATP Testing

Enable ATP products you would like to appear in the Device Type drop-down field when performing testing.

MicroSnap™

Toggle this on if you would like to run MicroSnap tests. Once turned on you'll be able to select products in the MicroSnap

product line.

Enzymes

Toggle this on if you would like to run Enzyme tests. Once turned on you'll be able to select Hygiena enzyme tests.

Note: Microsnap and Enzymes application will not appear unless enabled.

SureTrend Cloud

1. Login to [SureTrend Cloud](#).
2. Navigate to the **Instrument tab**.
3. Select the instrument you would like to manage products for.
4. Select **Settings** and then navigate to **Product Selections**.

The screenshot displays the SureTrend Cloud interface. At the top, the logo 'SureTrend by Hygiena' is visible. The navigation bar includes 'Dashboard', 'Results', 'Reports', and 'Instrument'. The 'Instrument' tab is active, showing a table of instruments. The selected instrument is 'Testing' with serial number 38019, site 'testing', and status 'Calibrated'. Below the instrument details, there are tabs for 'Locations/Samples', 'Plans', 'Users', 'Settings', and 'Calibration Checks'. The 'Settings' tab is selected, leading to the 'Product Selections' section. This section contains a table with columns for 'ATP', 'Lower', and 'Upper' values. The table lists four product types: UltraSnap, SuperSnap, AquaSnap Total, and AquaSnap Free, each with a checked checkbox and values of 20 for 'Lower' and 60 for 'Upper'. Below this, there are sections for 'Microorganisms By Sample' and 'Enzymes'.

SureTrend™ by Hygiena

Logout Settings Store Help

Dashboard Results Reports **Instrument**

Instruments

	Serial No	Instrument Name	Site	Calibration Check	Model	Software	Last Sync
	38019	Testing	testing	Calibrated	EnsureTouch	1.68.9	25-Sep-2020 10:47 AM

Instrument Name: Testing

Locations/Samples Plans Users **Settings** Calibration Checks

Product Selections

ATP	Lower	Upper
<input checked="" type="checkbox"/> UltraSnap	20	60
<input checked="" type="checkbox"/> SuperSnap	20	60
<input checked="" type="checkbox"/> AquaSnap Total	20	60
<input checked="" type="checkbox"/> AquaSnap Free	20	60

Microorganisms By Sample

Enzymes

5. Click the name of the product you would like to manage.

Product Selections

ATP

Microorganisms By Sample

Microorganisms enabled

MicroSnap Total

MicroSnap EB

MicroSnap Coliform

MicroSnap E. coli

Enzymes

6. If the product is disabled, check the box to enable it.

Product Selections

ATP

Microorganisms By Sample

Microorganisms enabled

MicroSnap Total

MicroSnap EB

MicroSnap Coliform

MicroSnap E. coli

Enzymes

7. Select the test device you would like to enable.

8. Sync the S/N you selected and the new product selections will be reflected on your instrument.

If you need assistance with product selection please contact Hygiena Technical Support by clicking [here](#).