



# How to Register Your EnSURE Touch

1. Log into [SureTrend Cloud](#).
2. Click the Register Instrument.


**Note:** If you do not see the Welcome page, select Settings > Welcome.

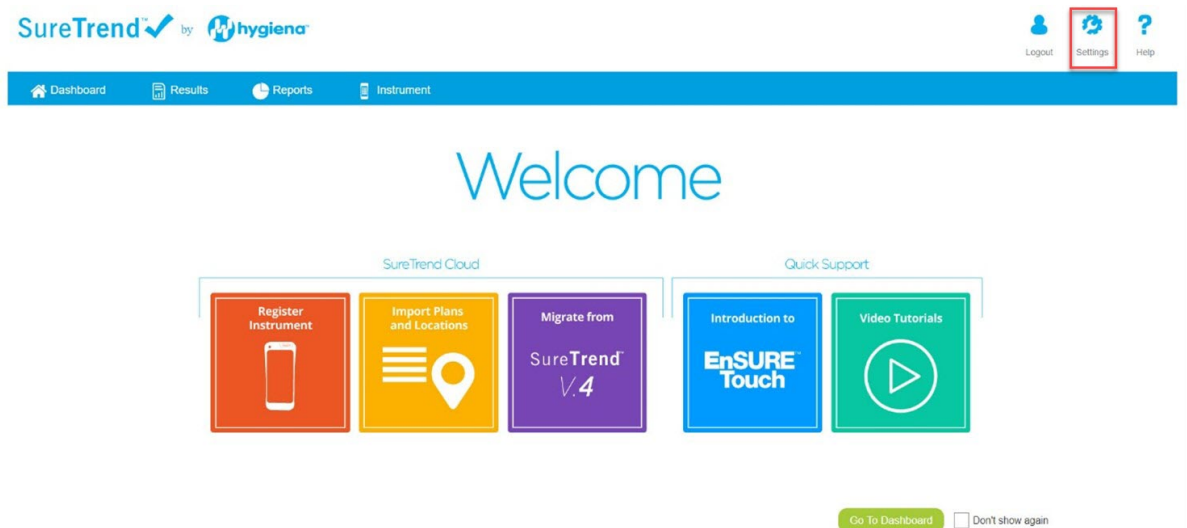


3. Name your EnSURE Touch and click Next. The name can be used to differentiate it from other instruments on the account. If you have multiple sites on your account, select the site to which your EnSURE Touch should be registered.
4. The cloud will display a registration token that you will use later in the registration process.

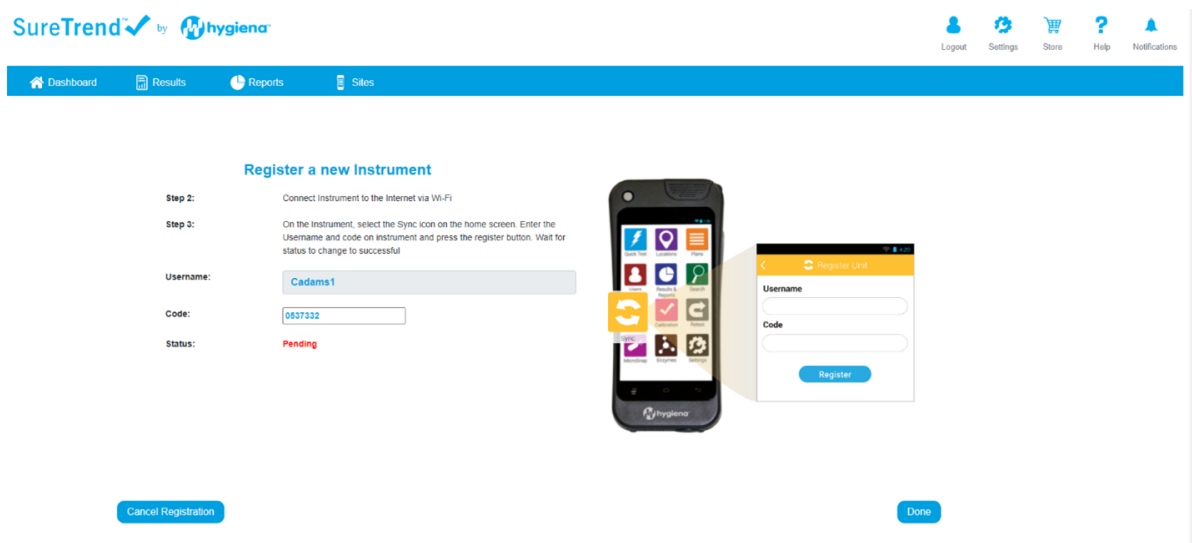

5. On your EnSURE Touch, open the Sync app.
6. Enter the Username and Code from SureTrend Cloud in the Sync app.

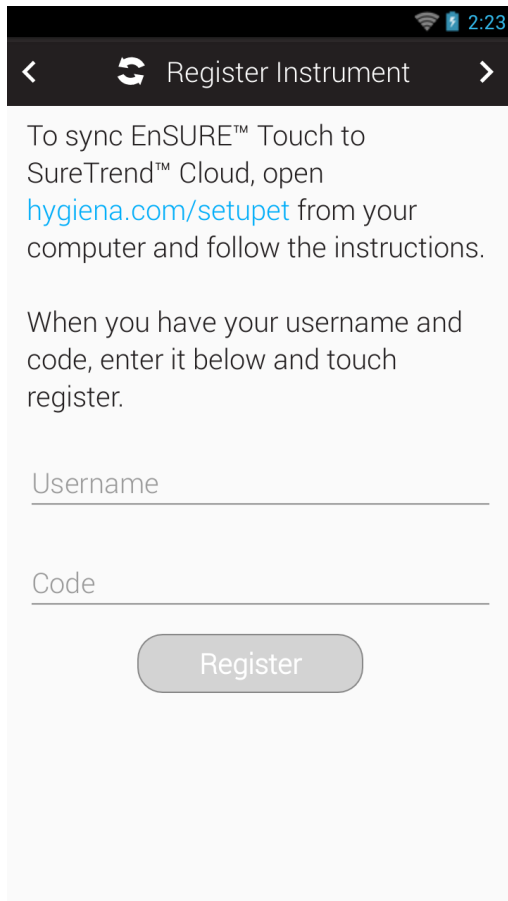

7. Once you have successfully Registered your Instrument your registration token in SureTrend Cloud will change from Pending to Successful, then select Done.

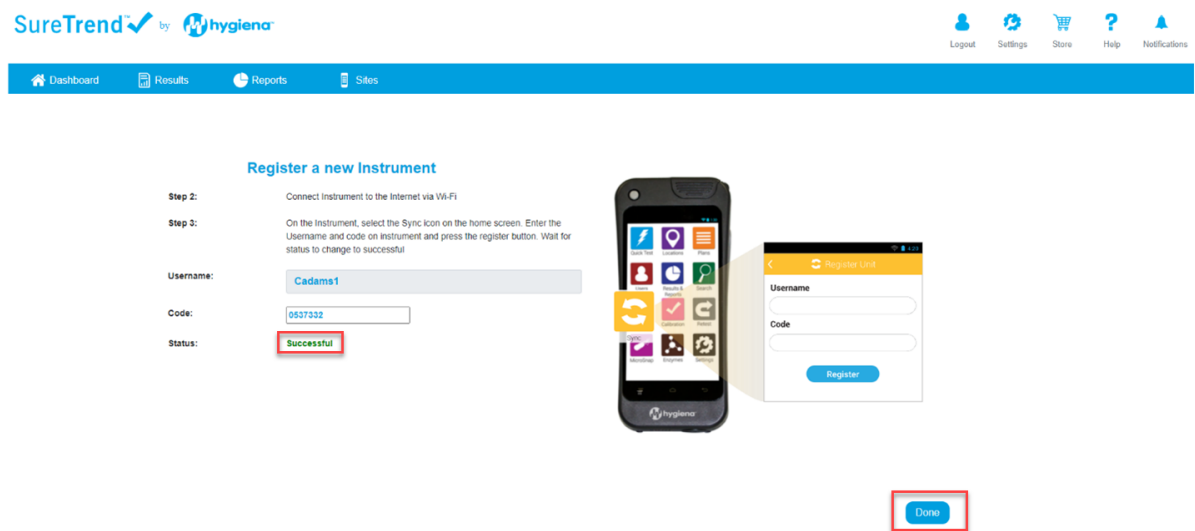

**Contact Hygiena Technical Support for further assistance.**

- Phone: 1-888-HYGIENA, option 2 (1-888-494-4362)
- Email: [techsupport@hygiena.com](mailto:techsupport@hygiena.com)
- [Submit a Support Ticket](#)

- [Schedule a Microsoft Teams meeting](#)