

# How to Register Your EnSURE Touch

## Registration

1. Log in to [SureTrend Cloud](#).
2. Click [here](#) to open to the registration page.
3. Enter a name your EnSURE Touch (optional).  
*Note: The name can be used to differentiate it from other instruments on the account.*
4. Select a site to register the EnSURE Touch if you have multiple sites.
5. Click Next.
6. On your EnSURE Touch, you will be prompted to register during the step up. Enter the Username and Code as it appears in SureTrend Cloud.  
*Note: If you have completed the step process without registering, you can open the Sync app to enter a username and code.*
7. Once you have successfully Registered your Instrument your registration token in SureTrend Cloud will change from Pending to Successful, then select Done.

## Contact Hygiena Technical Support for further assistance.

- Phone: 1-888-HYGIENA (1-888-494-4362, option 2)
- Email: [techsupport@hygiena.com](mailto:techsupport@hygiena.com)
- [Submit a Support Ticket](#)
- [Schedule a Microsoft Teams meeting with support](#)