


Adding a New/Replacement EnSURE Touch to a Site


Introduction

This article will show you how to add a new or replacement EnSURE Touch to your SureTrend Cloud account.

Solution

1. Login to [SureTrend Cloud](#).
2. Navigate to the Instrument Tab.
3. Select  located on the right hand side above the Instruments table.
4. Enter an instrument name (optional), and select a Site. Note: If you only have a default site you will not need to select this.

If this is a replacement instrument be sure to select the proper site so your plans and locations will sync.

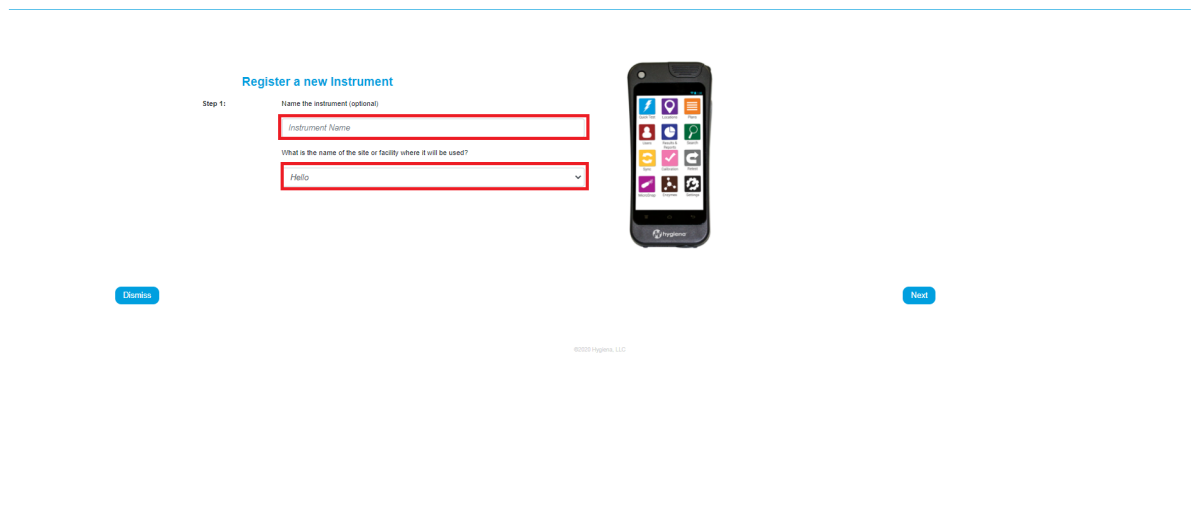

Register a new instrument

Step 1: Name the instrument (optional)

Instrument Name

What is the name of the site or facility where it will be used?

Hello

Dismiss Next

©2023 Vigorix, LLC

5. On the EnSURE Touch tap the Sync App and when prompted enter the Username and Code.

Register a new Instrument


Step 2: Connect Instrument to the Internet via Wi-Fi

Step 3: On the Instrument, select the Sync icon on the home screen. Enter the Username and code on instrument and press the register button. Wait for status to change to successful

Username:

Code:

Status: Pending

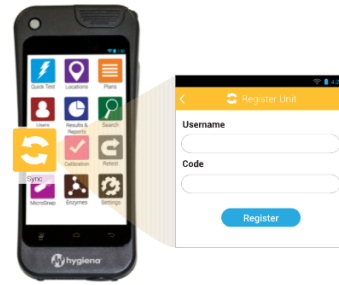

Cancel Registration

Done

6. Tap the sync now button to add your existing plans and locations.
Note: Plans and Locations are linked by the site the instrument are assigned to.
7. Your EnSURE Touch is now registered.