

# Adding Zone detail to Locations on the SureTrend Cloud and Ensure Touch

When creating a plan there is an option to add a Zone to each location. This will allow data in the Results tables to be viewed by Zone type.

The Zone is based on HACCP Zone criteria and as such there are 4 Zone options.


Zone 1 = Product Contact Surface

Zone 2 = Non PCS within close proximity to PCS

Zone 3 = Non PCS within process area but more removed from PCS

Zone 4 = Non PCS outside processing areas

On the Ensure Touch instrument, the Zone description is shown in the Zone text box.

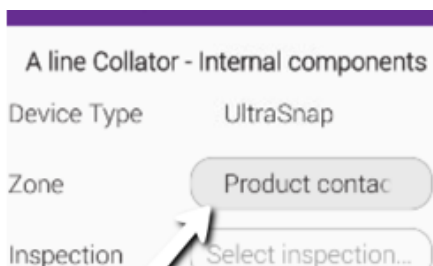


Zone detail can be added to locations either on the Sure Trend Cloud or on the Ensure Instrument.

## Enabling the Zone option

Before Zone detail can be added to locations the Zone feature needs to be enabled on the Sure Trend Cloud and Ensure Touch instrument.


## To enable the Zone feature on the Sure Trend Cloud.

Select the location tab and then settings. Scroll down to location settings and tick the enable zone box.

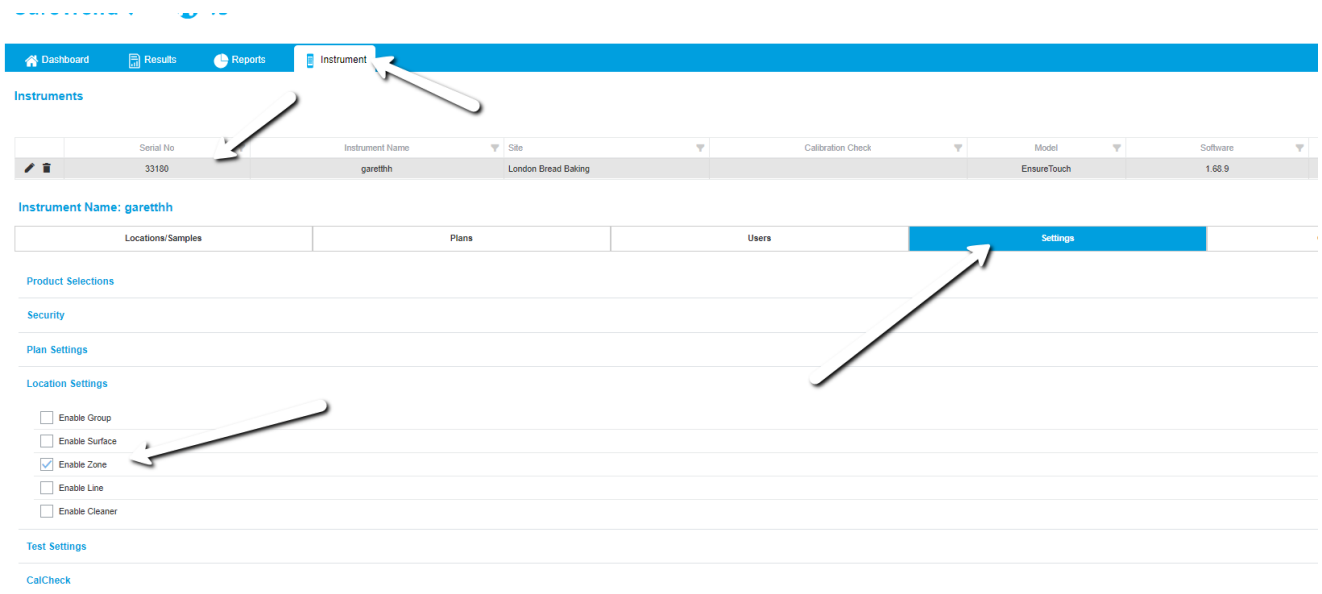


## To enable the Zone feature on the Ensure Touch

On the E Touch instrument from the man menu follow the steps below

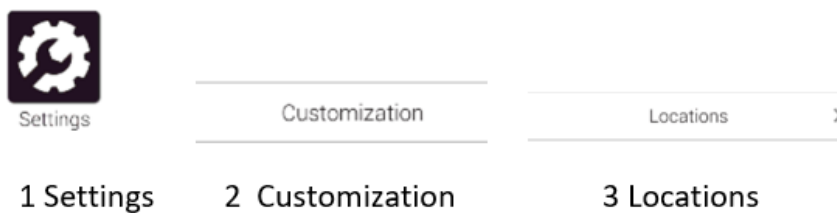

This will open the customization options. Then slide the Enable Zone button to on

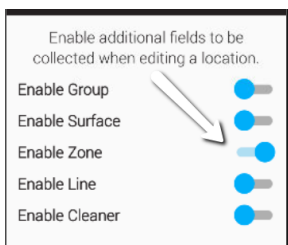


Once the Sure Trend Cloud and Ensure Touch instrument are enabled the Zones for each location can be added. Zones can be added to locations on the Sure Trend Cloud and Ensure Touch instrument.


## Setting Zones on Locations on the Sure Trend Cloud.

Select the Instrument Tab and select an instrument for the selected site.

Click the + icon.

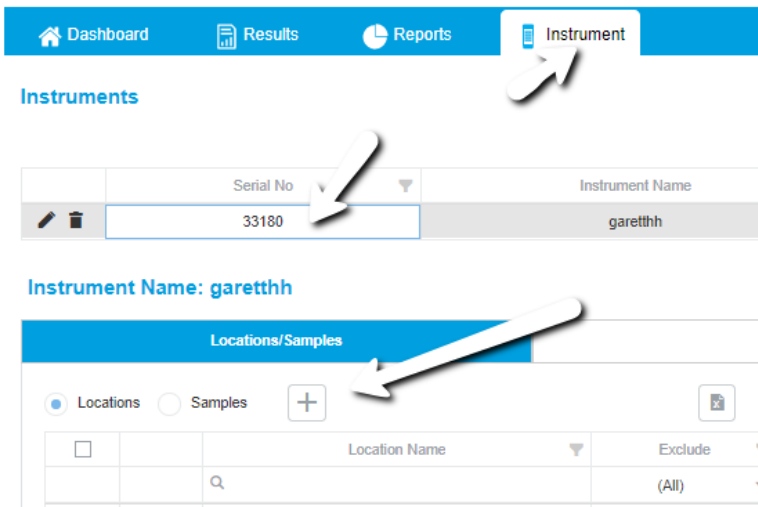

When Creating a new or Editing a location ( use the pencil icon to edit an existing location) add the Zone number to the Zone box

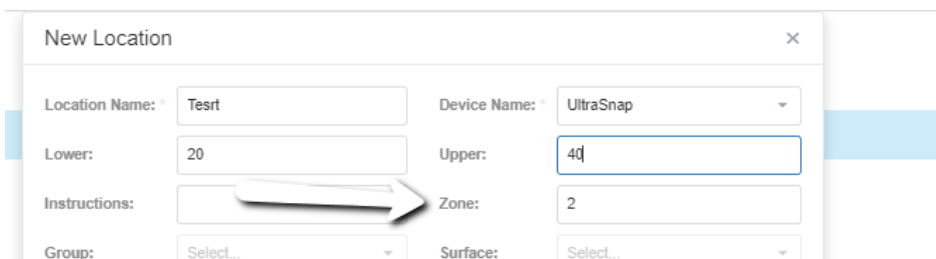

## Zone descriptions

**Note PCS = Product Contact Surface**

Zone 1 = Product Contact Surface

Zone 2 = Non PCS within close proximity to PCS


Zone 3 = Non PCS within process area but more removed from PCS

Zone 4 = Non PCS outside processing areas

Plan Name	Location	Device Name	Lower	Upper	Group	Surface	Zone
Zone Plan	Zone Test 1	UltraSnap	20	40			1
Zone Plan	Zone Test 2	UltraSnap	20	40			2
Zone Plan	Zone Test 3	UltraSnap	20	40			3
Zone Plan	Zone Test 4	UltraSnap	20	40			4
Zone Plan	Zone Test 5	UltraSnap	20	40			1
Zone Plan	Zone Test 6	UltraSnap	20	40			2
Zone Plan	Zone Test 7	UltraSnap	20	40			3
Zone Plan	Zone Test 8	UltraSnap	20	40			4
Zone Plan	Zone Test 9	UltraSnap	20	40			1
Zone Plan	Zone Test 10	UltraSnap	20	40			2


There are only 4 Zone types ( see above) so only numbers 1 to 4 can be added.

### Adding Zones on the E Touch.

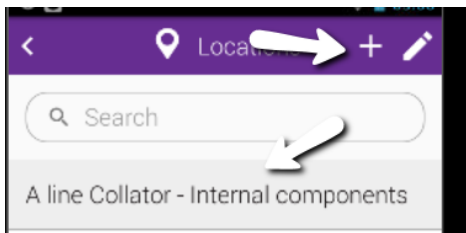


Select the Locations app on the Main menu-from here a new Location can be added + or an existing one edited using the pencil icon.

Tap on the Zone box to add the Zone criteria

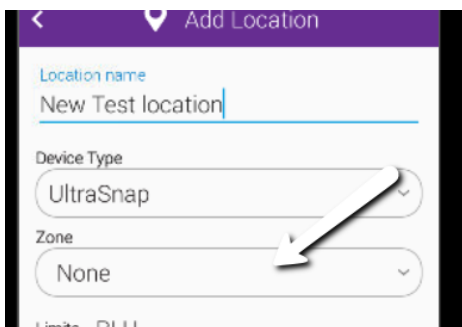


Select and tap the required Zone for the location. Note on the instrument the zones are not listed by number but the full HACCP description.

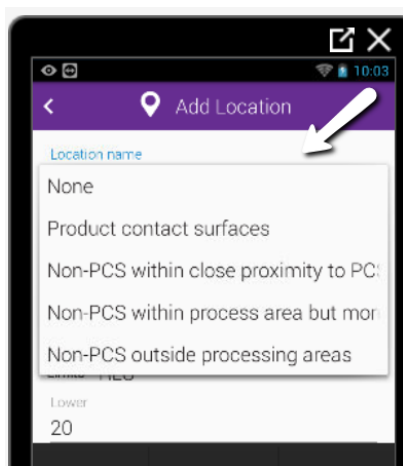

The Zone criteria will be added to the location - Select the back arrow to save.

When a location is then selected for testing the Zone criteria will appear and scroll across the zone box as shown below.

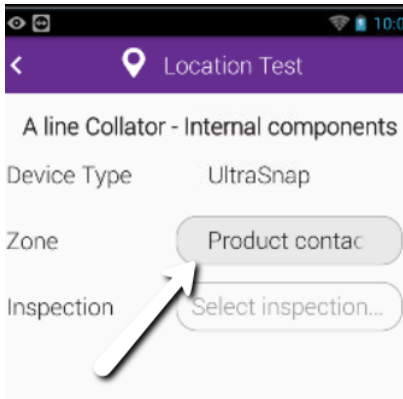

## Viewing results for Zones

Zone details will be found in the Results tables in the Zone Column. These can be filtered using the column filter.

The screenshot shows a web dashboard with a blue navigation bar containing 'Dashboard', 'Results', 'Reports', and 'Instrument'. Below the navigation bar, there are two dropdown menus: 'Type' set to 'ATP' and 'Period' set to 'This Year'. Below these is a text prompt: 'Drag a column header here to group by that column'. A table is displayed with the following data:

Location Name	Test Device	Lower	Upper	Notes	Zone	Rlu
Wine barrel	UltraSnap	20	40		Non-PCS within cl...	0

A white arrow points to the 'Zone' column header in the table.

The columns can then be filtered by the Zone type.

If you need any further support click on the icon below.



