

# Product Selection instructions for EnSURE Touch


## Introduction

Product Selections allows you to choose what apps and test devices are available on the EnSURE Touch™. In addition to testing ATP the EnSURE Touch also allows you test Micosnap and Enzymes. This article will show you how enable or disable products from EnSURE Touch and SureTrend Cloud.


## Solution

### EnSURE Touch


1. Touch the **Settings** App from the home screen. Next, touch **Customization**.

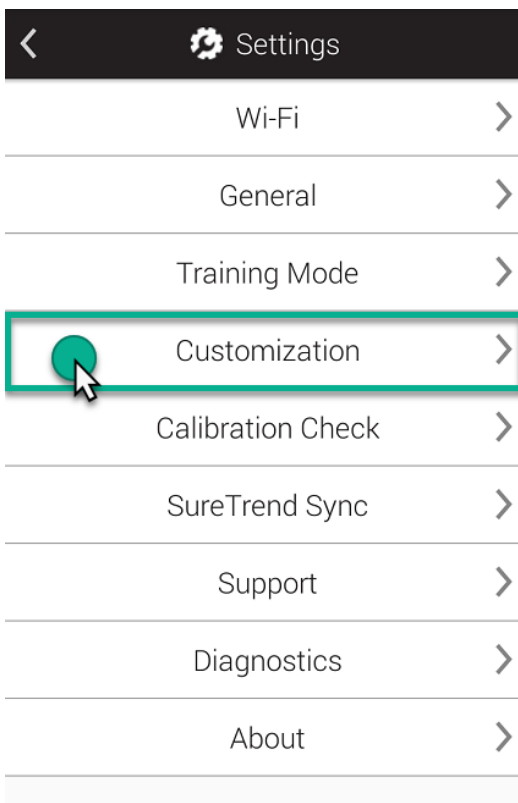

2. Select **Product Selections** □.

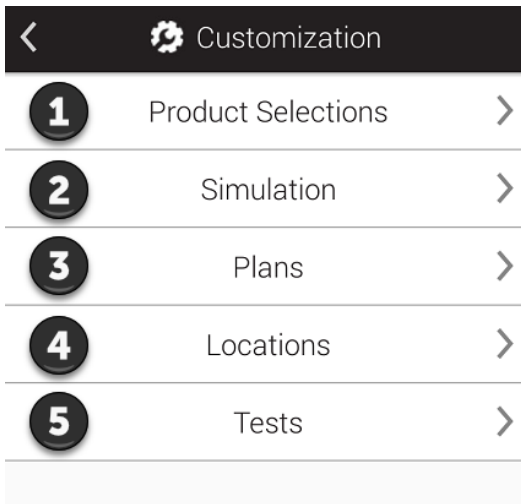

3. Select the product(s) you would like to enable/disable.

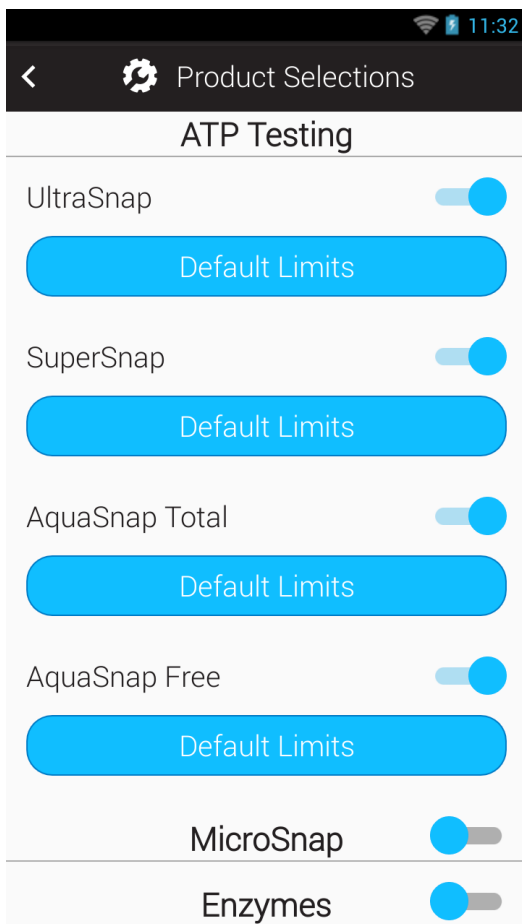

### **ATP Testing**

Enable ATP products you would like to appear in the Device Type drop-down field when performing testing.

### **MicroSnap™**

Toggle this on if you would like to run MicroSnap tests. Once turned on you'll be able to select products in the MicroSnap

product line.

## Enzymes

Toggle this on if you would like to run Enzyme tests. Once turned on you'll be able to select Hygiena enzyme tests.

**Note:** Microsnap and Enzymes application will not appear unless enabled.

# SureTrend Cloud

1. Login to [SureTrend Cloud](#).
2. Navigate to the **Instrument tab**.
3. Select the instrument you would like to manage products for.
4. Select **Settings** and then navigate to **Product Selections**.

The screenshot displays the SureTrend Cloud interface. At the top, the logo 'SureTrend by Hygiena' is visible on the left, and navigation links for Logout, Settings, Store, and Help are on the right. Below the logo is a blue navigation bar with tabs for Dashboard, Results, Reports, and Instrument. The 'Instrument' tab is selected. Underneath, the 'Instruments' section shows a table with one instrument: Serial No. 38019, Instrument Name: Testing, Site: testing, Calibration Check: Calibrated, Model: EnsureTouch, Software: 1.68.9, and Last Sync: 25-Sep-2020 10:47 AM. Below the table, the 'Instrument Name: Testing' section has a sub-navigation bar with tabs for Locations/Samples, Plans, Users, Settings (selected), and Calibration Checks. The 'Product Selections' section contains a table with columns for ATP, Lower, and Upper. The table lists four products: UltraSnap, SuperSnap, AquaSnap Total, and AquaSnap Free, each with a checked checkbox and values of 20 for Lower and 60 for Upper. At the bottom, there are sections for 'Microorganisms By Sample' and 'Enzymes'.

**SureTrend™** by **Hygiena**

Logout Settings Store Help

Dashboard Results Reports Instrument

Instruments

|  | Serial No | Instrument Name | Site    | Calibration Check | Model       | Software | Last Sync            |
|--|-----------|-----------------|---------|-------------------|-------------|----------|----------------------|
|  | 38019     | Testing         | testing | Calibrated        | EnsureTouch | 1.68.9   | 25-Sep-2020 10:47 AM |

Instrument Name: Testing

Locations/Samples Plans Users Settings Calibration Checks

Product Selections

| ATP  | Lower | Upper |
|--|-------|-------|
| <input checked="" type="checkbox"/> UltraSnap      | 20    | 60    |
| <input checked="" type="checkbox"/> SuperSnap      | 20    | 60    |
| <input checked="" type="checkbox"/> AquaSnap Total | 20    | 60    |
| <input checked="" type="checkbox"/> AquaSnap Free  | 20    | 60    |

Microorganisms By Sample

Enzymes

5. Click the name of the product you would like to manage.

#### Product Selections

ATP

**Microorganisms By Sample**

Microorganisms enabled

MicroSnap Total

MicroSnap EB

MicroSnap Coliform

MicroSnap E. coli

Enzymes

6. If the product is disabled, check the box to enable it.

#### Product Selections

ATP

**Microorganisms By Sample**

Microorganisms enabled

MicroSnap Total

MicroSnap EB

MicroSnap Coliform

MicroSnap E. coli

Enzymes

7. Select the test device you would like to enable.

8. Sync the S/N you selected and the new product selections will be reflected on your instrument.

If you need assistance with product selection please contact Hygiena Technical Support by clicking [here](#).