


Modifying a location


Step One

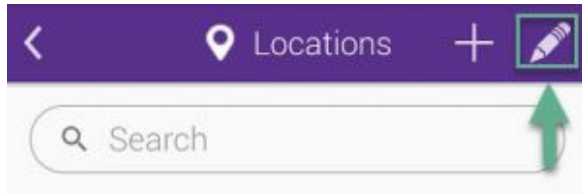

[Click to enlarge](#)

Touch on the pencil icon as shown in the image to modify an existing location.

Step Two

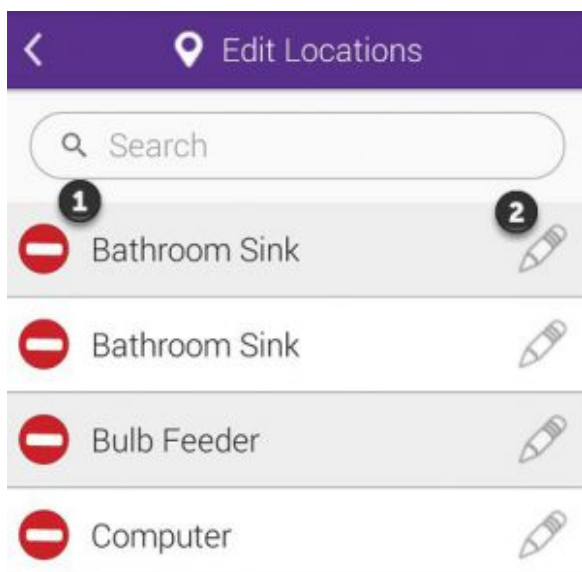


[Click to enlarge](#)

Touch the pencil icon as shown at

2

to begin modifying the location.

Step Three

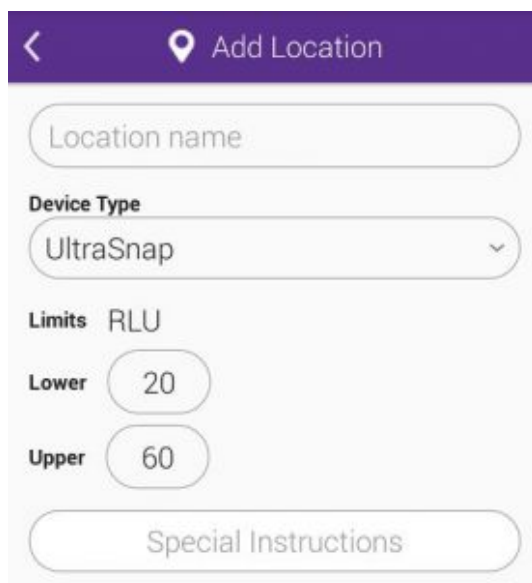

The screenshot shows a mobile application interface for adding a location. At the top is a purple header bar with a back arrow on the left and a location pin icon followed by the text 'Add Location'. Below the header is a form with several fields: a text input field labeled 'Location name'; a dropdown menu labeled 'Device Type' with 'UltraSnap' selected; a section titled 'Limits' containing two sub-sections, 'RLU' and 'RLU', each with a 'Lower' and 'Upper' value (20 and 60 respectively); and a text input field at the bottom labeled 'Special Instructions'.

[Click to enlarge](#)

You will be able to change the location's name, device type and RLU limits. You'll also be able to add or change any special instructions.