


# Modifying a location


## Step One

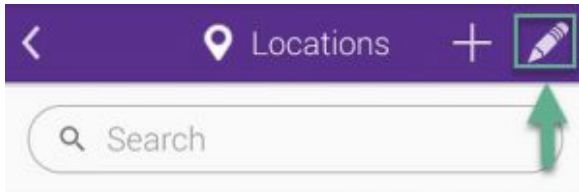

Click to enlarge

Touch on the pencil icon as shown in the image to modify an existing location.

## Step Two

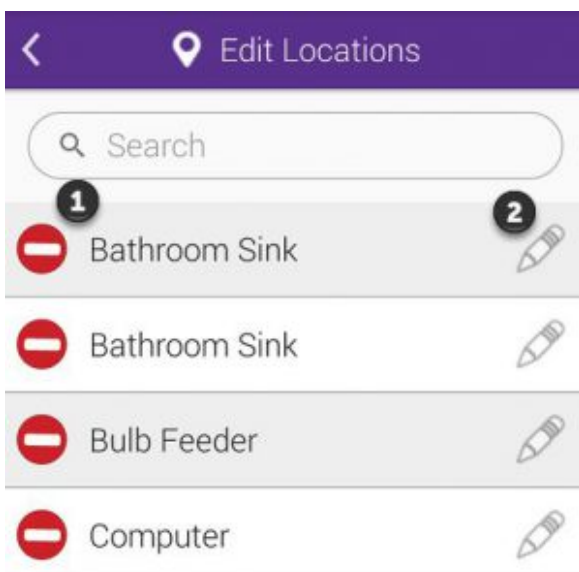


Click to enlarge

Touch the pencil icon as shown at

2

to begin modifying the location.

## Step Three

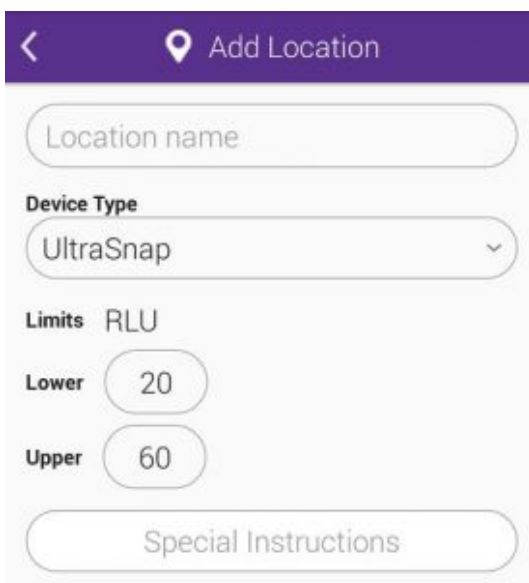

The screenshot shows a mobile application interface for adding a location. At the top, there is a purple header with a back arrow on the left and a location pin icon followed by the text "Add Location". Below the header is a form with several fields: a text input field labeled "Location name", a dropdown menu labeled "Device Type" with "UltraSnap" selected, a section titled "Limits" with "RLU" as a sub-label, two numeric input fields labeled "Lower" (with the value 20) and "Upper" (with the value 60), and a text input field at the bottom labeled "Special Instructions".

[Click to enlarge](#)

You will be able to change the location's name, device type and RLU limits. You'll also be able to add or change any special instructions.