


Modifying a location


Step One

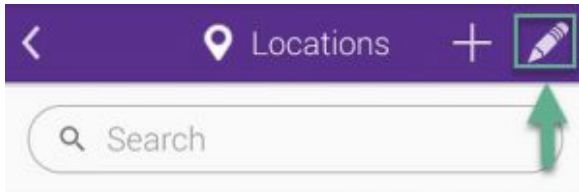

Click to enlarge

Touch on the pencil icon as shown in the image to modify an existing location.

Step Two

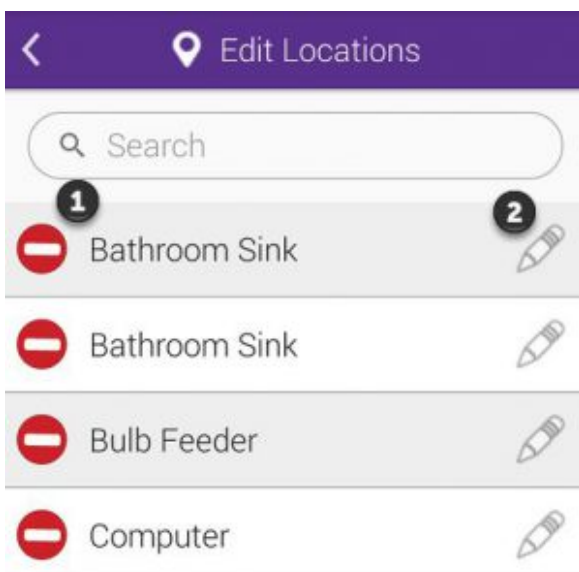


Click to enlarge

Touch the pencil icon as shown at

2

to begin modifying the location.

Step Three

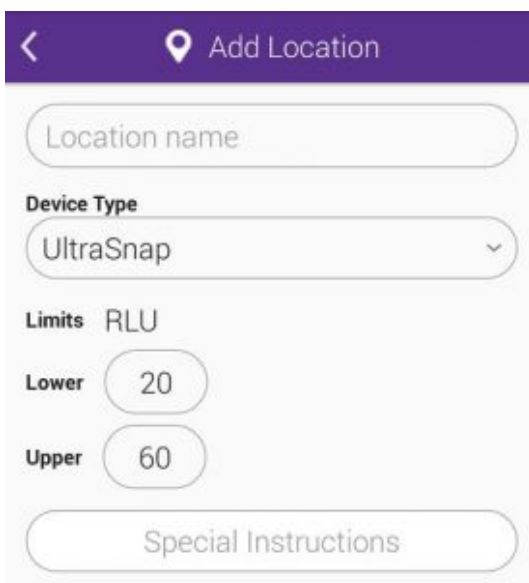

The screenshot shows a mobile application interface for adding a location. At the top, there is a purple header with a back arrow on the left and a location pin icon followed by the text "Add Location". Below the header is a form with several fields: a text input field for "Location name", a dropdown menu for "Device Type" currently showing "UltraSnap", a section for "Limits RLU" with two numeric input fields: "Lower" set to 20 and "Upper" set to 60, and a text input field for "Special Instructions".

[Click to enlarge](#)

You will be able to change the location's name, device type and RLU limits. You'll also be able to add or change any special instructions.