



# Running tests from a plan

## Step One

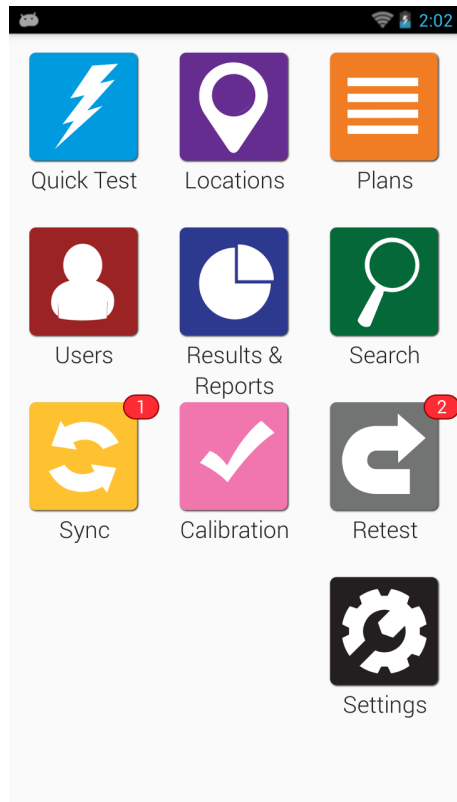

From the **Home Screen**, tap the **Plans** app.

# Step Two

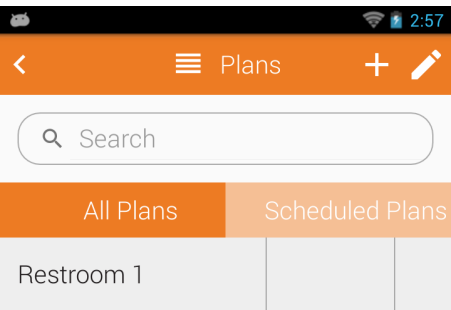

Tap the name of the plan that you want to test.

# Step Three

2:57

< Plan Locations >

Restroom 1

Started 08/26/2020 at 2:57 PM

Plan Locations Tested 0 of 3

Restroom Door Handle	
Restroom Sink	
Restroom Stall Door	

To run the tests in order, tap the **Start** button.

## Step Four

Restroom 1

Restroom Door Handle

Location 1 of 3

Device Type UltraSnap

Group

Surface

Zone None

Line

Cleaner

Run Test

Insert your test device and tap the **Run Test** button.

## Step Five

After the test is finished, tap the **Next** button to proceed to the next test.

## Step Six

Repeat Steps 4 & 5 for each location in the plan. Tap the **Done** button when the last test is finished.