


# Chapter 3 | Results

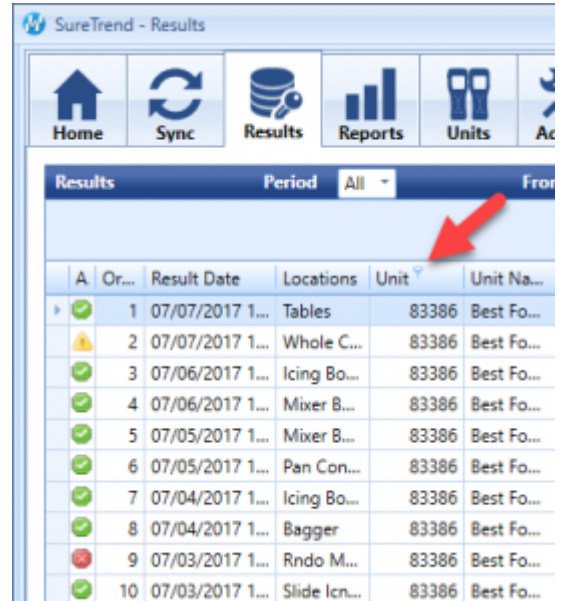

The screenshot shows the 'SureTrend - Results' interface. At the top, there is a navigation bar with icons for Home, Sync, Results, Reports, and Units. Below this is a 'Results' section with a 'Period' dropdown set to 'All' and a 'From' field. A table of results is displayed below, with columns: A, Or..., Result Date, Locations, Unit, and Unit Na... A red arrow points to the small pin icon next to the 'Unit' column header. The table contains 10 rows of data, each with a status icon (green checkmark or yellow warning triangle), a number, a date, a location, a unit number, and a unit name.

| A | Or... | Result Date | Locations       | Unit         | Unit Na...       |
|---|-------|-------------|-----------------|--------------|------------------|
| ▶ | ✓     | 1           | 07/07/2017 1... | Tables       | 83386 Best Fo... |
|   | ⚠     | 2           | 07/07/2017 1... | Whole C...   | 83386 Best Fo... |
|   | ✓     | 3           | 07/06/2017 1... | Icing Bo...  | 83386 Best Fo... |
|   | ✓     | 4           | 07/06/2017 1... | Mixer B...   | 83386 Best Fo... |
|   | ✓     | 5           | 07/05/2017 1... | Mixer B...   | 83386 Best Fo... |
|   | ✓     | 6           | 07/05/2017 1... | Pan Con...   | 83386 Best Fo... |
|   | ✓     | 7           | 07/04/2017 1... | Icing Bo...  | 83386 Best Fo... |
|   | ✓     | 8           | 07/04/2017 1... | Bagger       | 83386 Best Fo... |
|   | ✗     | 9           | 07/03/2017 1... | Rndo M...    | 83386 Best Fo... |
|   | ✓     | 10          | 07/03/2017 1... | Slide Icn... | 83386 Best Fo... |

## Filtering Columns


Clicking on the little pin next to any column title will allow you to filter that column by any data present in that column.

With this feature you can customize the results view to only show the exact data you are interested in, i.e. showing results from specific units, users, locations or dates only.

## View custom time periods

You can select to view only results from a specific period here. For example: this month, last month, this week or last week. You can choose All to display all results and reset any custom time periods you may have turned on.

You can also use the From and To to show custom date ranges of your choosing. Just select the starting and ending dates and view all results within that time frame.

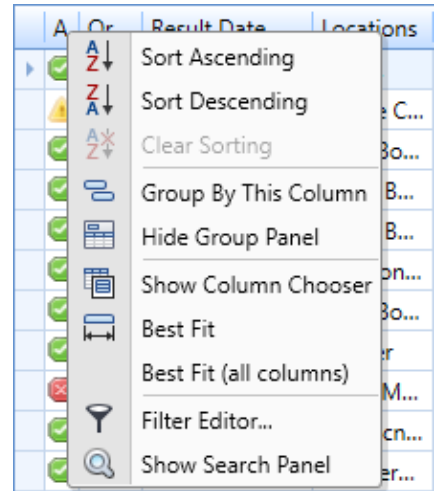

Right click on any column title and to choose from a menu of options that will allow you to further customize the view of your results. You can even choose to resize columns to make them easier to view.