


Chapter 4 | Reporting

Report Settings

- [Filter](#)
- [Specify](#)
- [Save](#)
- [Report Properties](#)
- [Display Options](#)
- [Color Options](#)

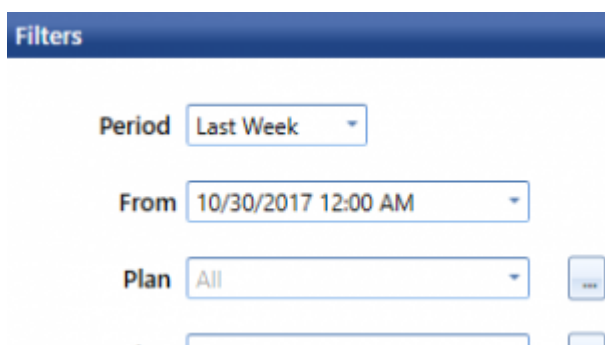


The screenshot shows a 'Filters' section with the following controls:


- Period:** A dropdown menu currently set to 'Last Week'.
- From:** A date and time selector currently set to '10/30/2017 12:00 AM'.
- Plan:** A dropdown menu currently set to 'All', with a small '...' button to its right.
- Below the 'Plan' dropdown, there is a partially visible checkbox and another dropdown menu.

Here you can choose what will be in your report. By default All is chosen which means all plans, all users, all locations, etc. will be included in the reports data.

Checking the box **Include Deleted Rows** will include deleted results in the report.

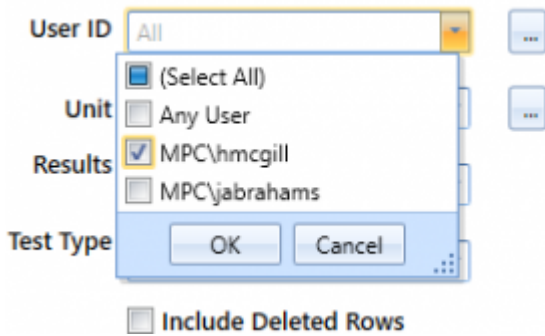


Clicking on any of the fields will allow you to pick and choose what data will appear in your reports. For example, you can specify specific plans, locations, users, or even units to include in your reports if you do not want to include all.

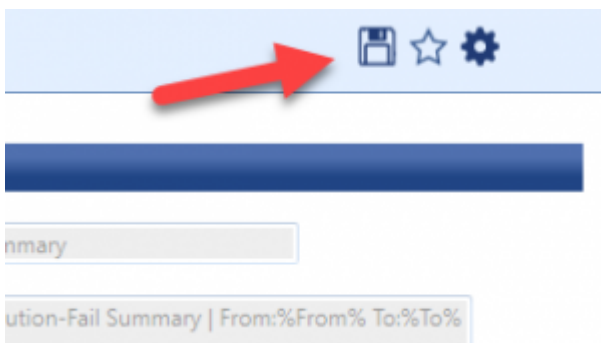

You can save your entered information if this is a report you conduct often by clicking the floppy disc icon.

Pressing the star will allow you to add this report to your Favorites list. Clicking the gear icon will allow you to save and customize this report.

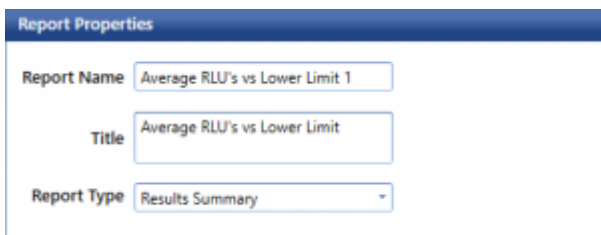

This panel is only available when working with a saved or custom report.

You can change the report name that appears in the Customized Reports list and

choose a Title that will appear at the top of the report.



Display Options


Chart Type: Bar




X Axis:

Y Axis:

This panel is only available when working with a saved or custom report.

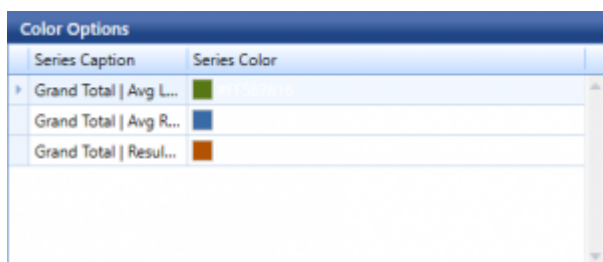Select the type of chart you want your report to include. You can select from a variety of bar, line, point and pie charts. You may also edit the axis text.



Series Caption	Series Color
Grand Total Avg L...	
Grand Total Avg R...	
Grand Total Resul...	

This panel is only available when working with a saved or custom report.

Select the color(s) you want to appear in your report charts, graphs, or tables.