


Product Selections

Customizations are optional. These settings may be changed any time without risk to your data.


Step One

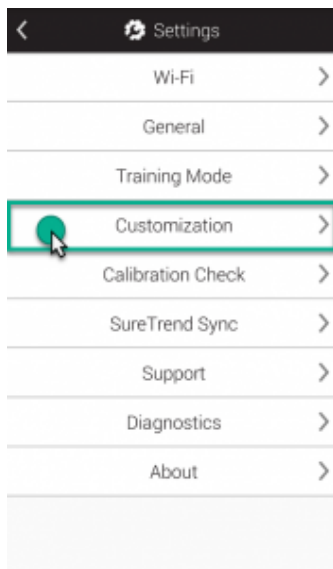

Click to enlarge

Touch on the **Settings** button from the Home screen. Next, touch **Customization**.


Step Two

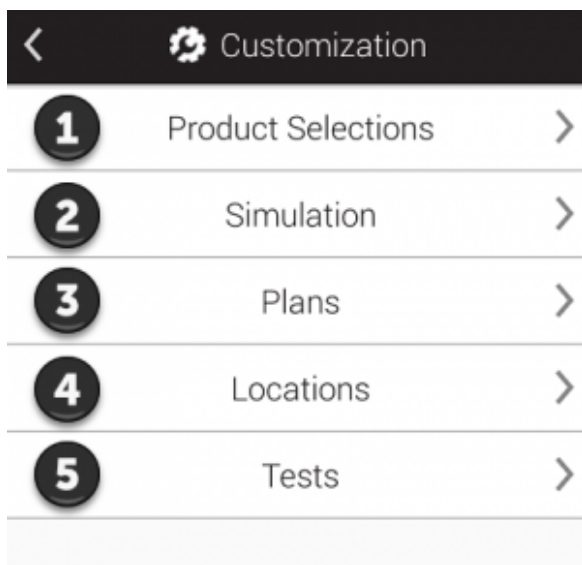

Click to enlarge

Touch **Product Selections** (as shown at **1** in the image).

Step Three

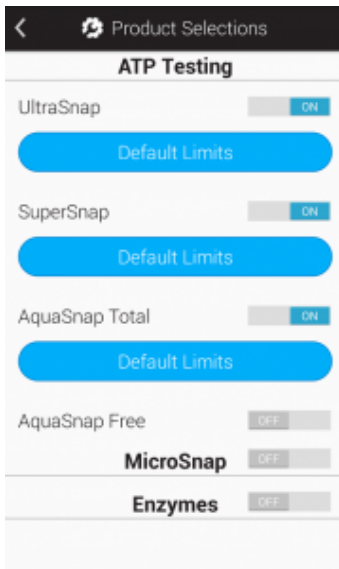

Click to enlarge

ATP Testing

Toggle on any products you would like to appear in the Device Type drop-down field when conducting regular tests and quick tests.

MicroSnap

Toggle this on if you'd like to run MicroSnap tests. Once toggled on you'll be able to select any specific surfaces or products in the MicroSnap line you want to use.

Enzymes

Toggle this on if you'd like to run Enzyme tests. Once toggled on you'll be able to select any specific tests you want to use.