

Chapter 5 | Units

Adding & Managing Units

There are two ways you can add a Unit to SureTrend.

1. Connect your unit to the machine and perform a sync. This will automatically import the unit and sync any data.
2. Click the **Plus** icon in the Units section (as shown below). In the window that pops up click **Discover** and then click **Add**. Be sure **Auto detect USB cable** is checked.




Click to enlarge



Click to enlarge

- 1
- 2
- 3
- 4
- 5

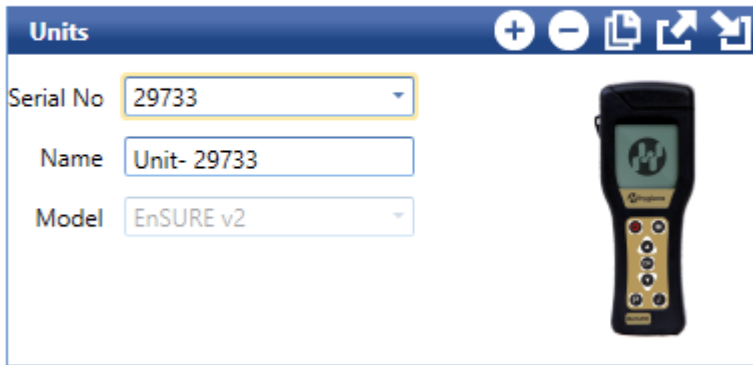

1	Add/Discover Unit: Add a non-existing unit to SureTrend
2	Delete Unit: Remove the selected unit from SureTrend
3	Copy Unit: Copy program settings from an existing unit to another unit
4	Export unit to XML file: Export a unit from SureTrend to an XML file
5	Import unit to XML file: Import a unit into SureTrend from an XML file