

# Adding & Managing Locations

Or...	Location Name	Gro...	Surf...	Up...	Lo...	R...	Zone	Room Nu...	Perso...	No...	Corrective A...
*	Grid New Row Text										
0	90 degree Belt			60	59	0	Product contact su...	0	0		
1	Accumulator Infeed			60	59	0	Product contact su...	0	0		
2	Accumulator Outf...			60	59	0	Product contact su...	0	0		
3	Any Location			60	59	0	Product contact su...	0	0		
4	Batter Holding Tank			60	59	0	Product contact su...	0	0		
5	Batter Mixing Tank			60	59	0	Product contact su...	0	0		

1	Select from your collection of locations
2	Paste locations from your clipboard
3	Add a new location
4	Remove the selected location(s)
5	Save and commit all changes
6	Quickly add new locations by typing their name here and pressing Enter when done
7	Click the pin icon in any column to filter what you wish to see

## How locations can be used

Usually locations are used as the physical locations you test with a unit. And since SureTrend can hold up to 5,000 locations you can add plenty and group them by using the Groups, Surfaces, or Room Number fields. You may also create custom

fields as described in the *Location Configuration* section of Chapter 6.