

# Transferring Unit Data

One of the ways you can transfer data is by copying it from a unit already in SureTrend. With the Copy Unit feature you can transfer program data (including users, locations, and plans) to another unit already in SureTrend or to a unit connected to SureTrend via USB cable.

Click **Discover** to transfer data to a unit connected via USB cable. To transfer to a unit already in SureTrend select it from the **Serial No.** drop-down box. Click **Copy** once you are ready to proceed and copy the data.

**1**

**Units**

Serial No. 29733

Name Unit- 29733

Model EnSURE v2



**2**

**From Unit**

Serial No. 29773

Name Unit- 29773

Model EnSURE v2



**To Unit**

☒ Auto detect USB cable

Serial No. 111111

Name Unit - 90036

Model SystemSURE Plus v2

Serial Port COM3

Discover

Copy Cancel

Once the copy is completed do a sync to transfer the data onto your unit.