


# Finding the MAC address


## Why would I need the MAC address?

Sometimes only specific, known devices are allowed onto a network. Other times a specific device needs to be identified on a network. When these situations arise you may need your EnSURE Touch's MAC address which is a string of characters that are unique and identifies it. DHCP servers use the MAC address to identify devices and give some fixed IP addresses. If your network environment requires needing your devices MAC address follow these instructions to find it.

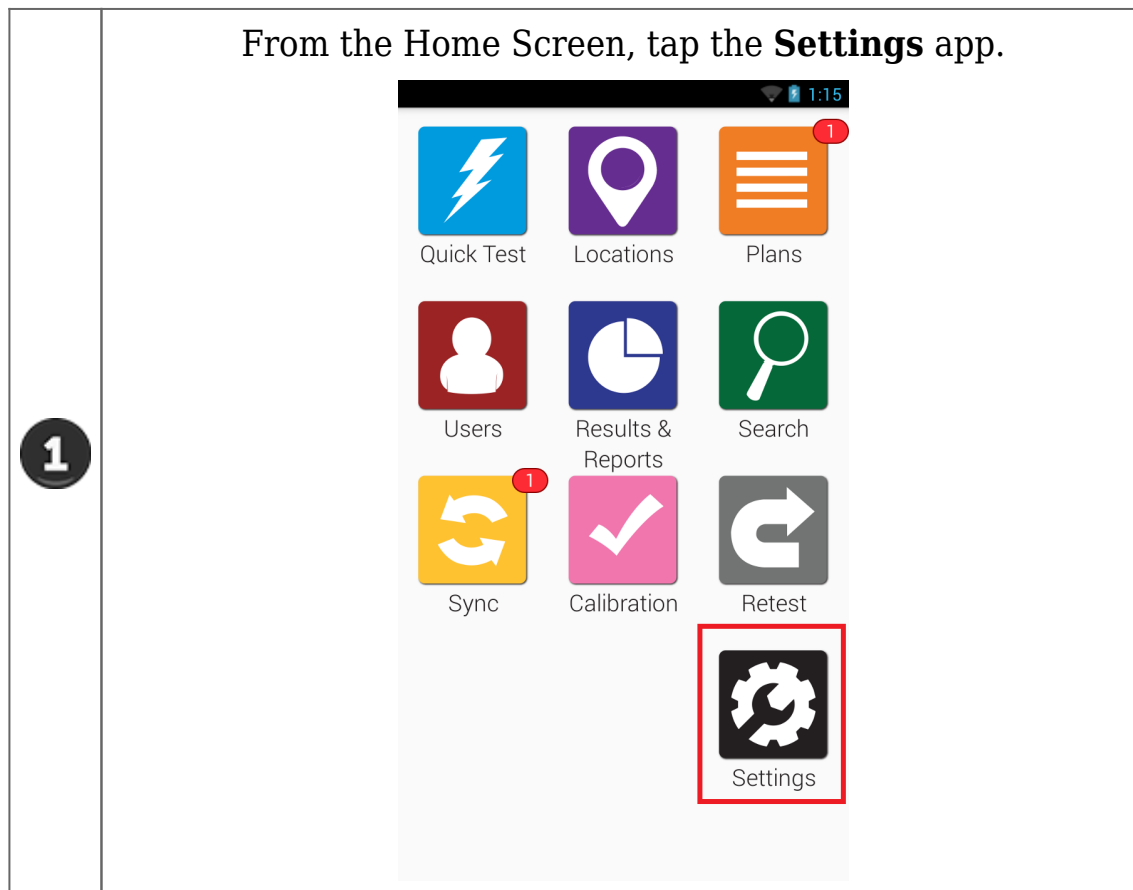

## Tap **About**.

2

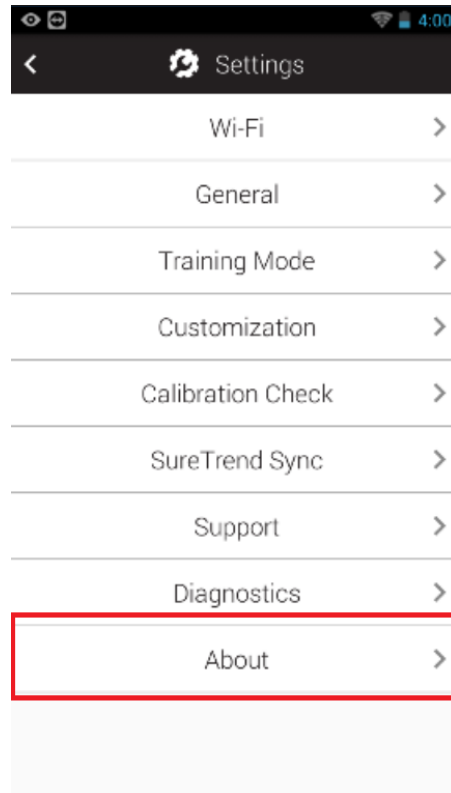

## Tap **Hardware**.

The instrument's MAC Address is the seventh item on the list.

3

