



Recording an Allergen Quick Test

Recording an Allergen Quick Test

1. Open the Allergen app.

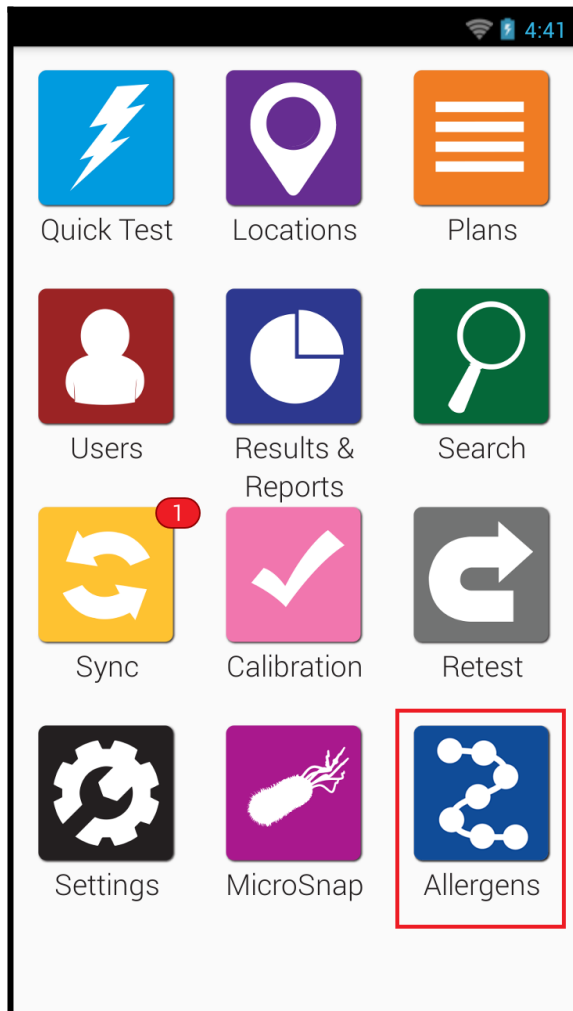

2. Tap Quick Test.

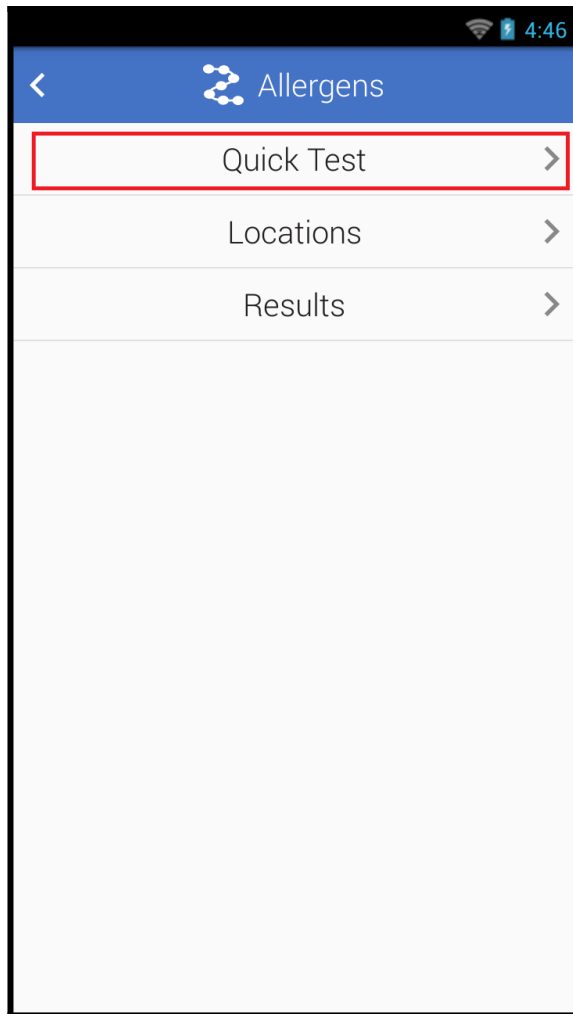







3. Select the correct test device from the Device Type dropdown menu.
4. For test devices with multiple allergen tests, select the specific allergen from the Allergens dropdown menu.

Quick Test

Device Type AlerTox Sticks

Allergens Total Milk

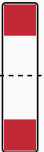Select the result that matched your Device

	Positive	Negative	Invalid
Control			
Positive			







5. Tap the correct result that matches your test result.

Quick Test

Device Type AlerTox Sticks

Allergens Total Milk

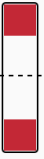Select the result that matched your Device

	Positive	Negative	Invalid
Control			
Positive			

Save

6. Tap the Save button.
7. Review the test information and tap the Done button to finish recording the result.

10:08

<

Quick Test Result

Device Type

AlerTox Sticks

Allergens

Total Milk

Time

10/20/2021 10:08 AM

Positive

Negative

Invalid

Control

Positive

Done