



Recording an Allergen Quick Test

Recording an Allergen Quick Test


1. Open the Allergen app.

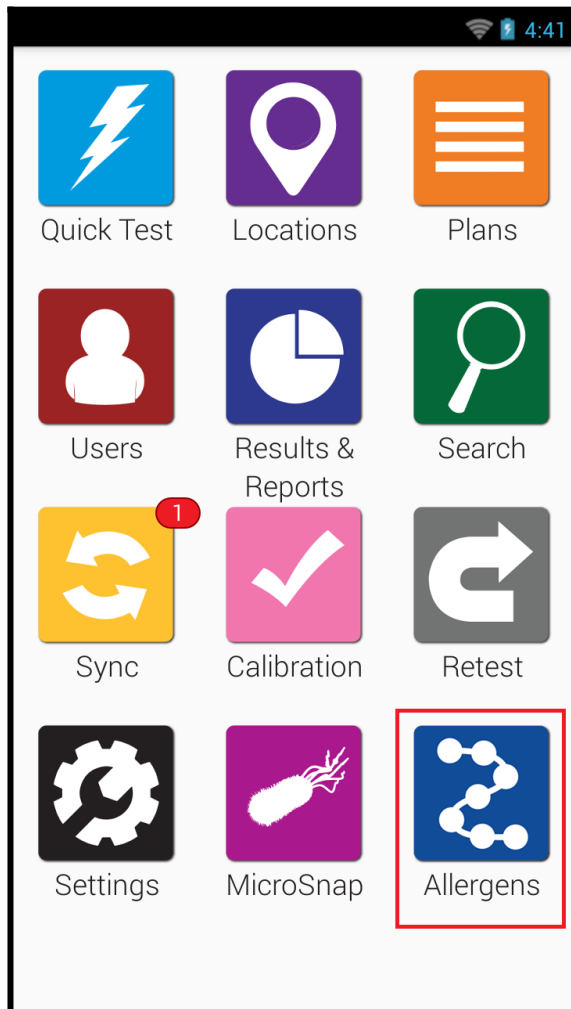


2. Tap Quick Test.

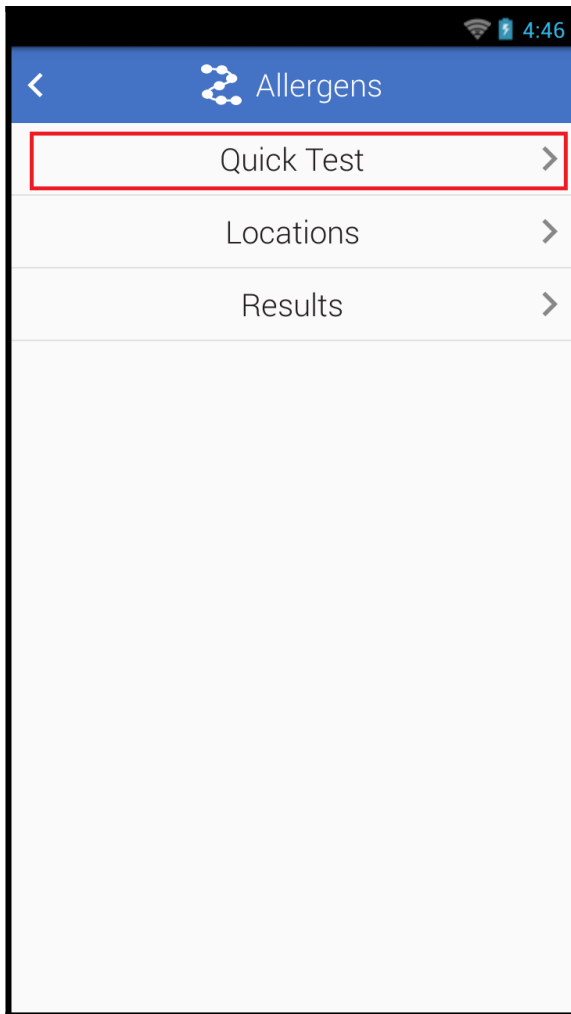

3. Select the correct test device from the Device Type dropdown menu.
4. For test devices with multiple allergen tests, select the specific allergen from the Allergens dropdown menu.

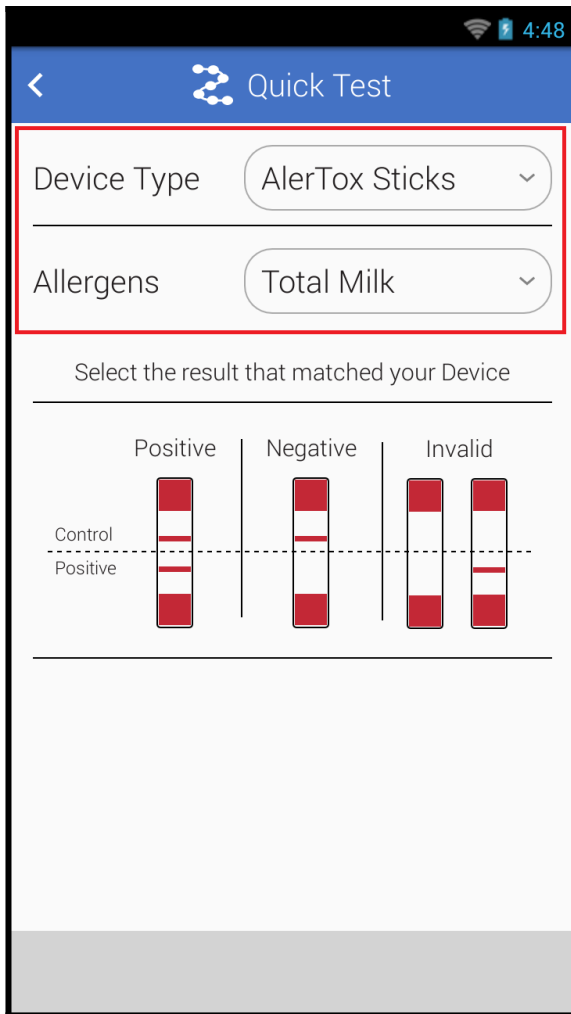

5. Tap the correct result that matches your test result.

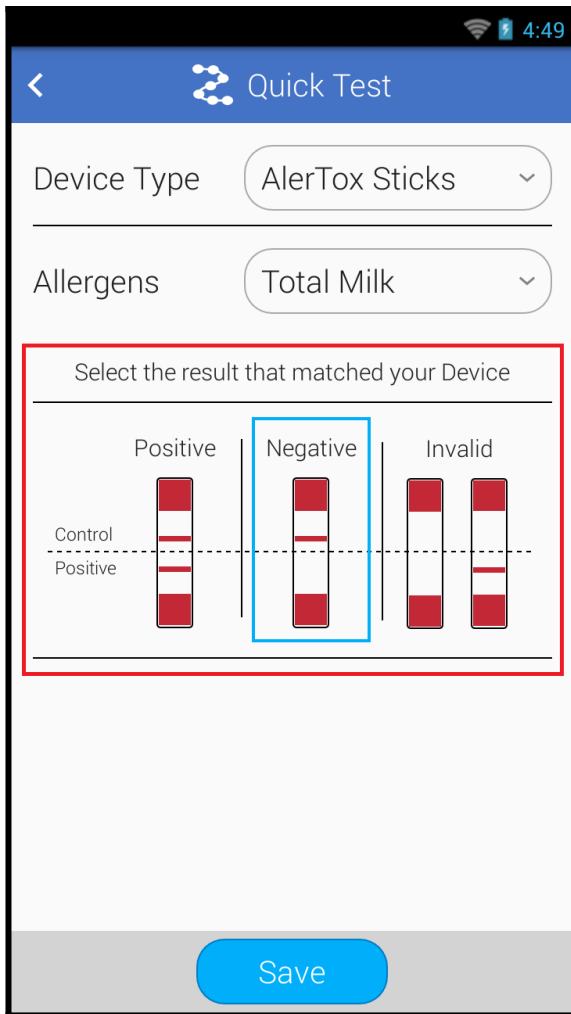


6. Tap the Save button.
7. Review the test information and tap the Done button to finish recording the result.

Quick Test Result

Device Type AlerTox Sticks

Allergens Total Milk

Time 10/20/2021 10:08 AM

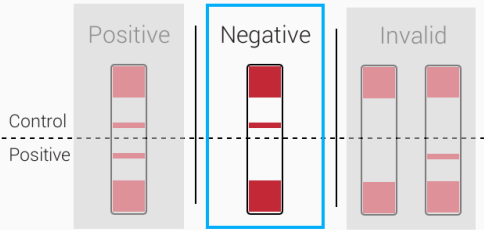

Done